



# Afla CLEAN™ SMART Handling

## 1. Sample

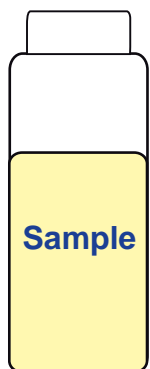


## 2. Apply Sample

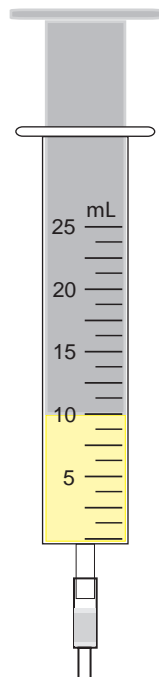

10 mL sample  
with 3 mL/min  
3,5 min


## 3. Wash Column

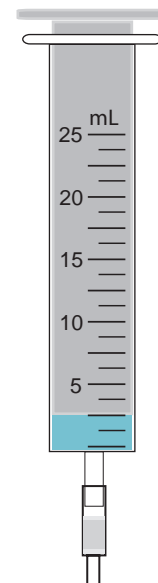

2 mL water  
with 3 mL/min  
0,5 min

→ Drying

## 4. Eluting Sample

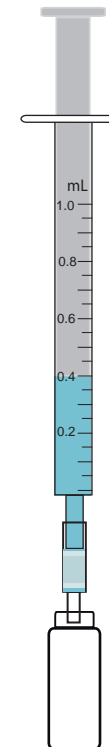


0,4 mL MeOH  
5 min waiting time


**IAC-Clean-up time < 9 min**

# Ota CLEAN™ SMART Handling

## 1. Sample

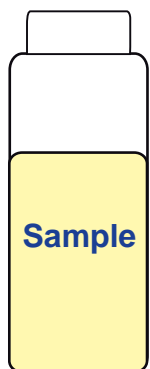


## 2. Apply Sample

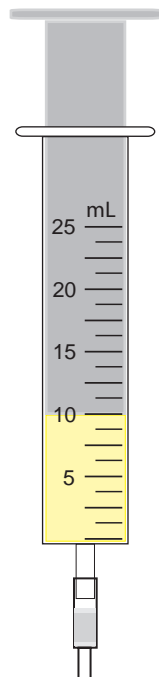

10 mL sample  
with 1,5 mL/min  
6,5 min


## 3. Wash Column

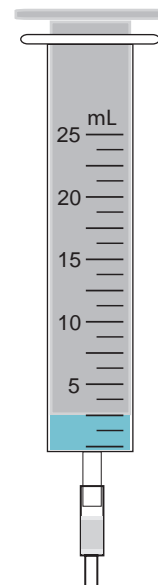

2 mL water  
with 1,5 mL/min  
1,5 min

→ Drying

## 4. Eluting Sample

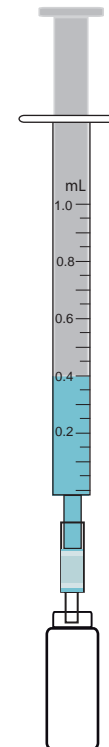

0,4 mL MeOH  
5 min waiting time

**IAC-Clean-up time < 13 min**