



Agilent Genomic DNA ScreenTape Assay Quick Guide for 4200 TapeStation System

The Agilent 4200 TapeStation system (G2991AA) is an automated platform for scalable, flexible, faster and more reliable electrophoresis. The Genomic DNA ScreenTape assay is designed for analyzing genomic DNA in the sizing range from 200 to >60000 bp.

Specifications

Analytical Specifications	Genomic DNA ScreenTape Assay
Sizing Range	200 to >60000 bp
Sensitivity ¹	0.5 ng/ μ L
Sizing Precision ²	200 – 15000 bp: 15 % CV
Sizing Accuracy ²	200 – 15000 bp: \pm 15 %
Quantitative Precision ³	15 % CV
Quantitative Accuracy ³	\pm 20 %
Quantitative Range	10 – 100 ng/ μ L
DIN functional range ⁴	5 – 300 ng/ μ L
Physical Specifications	
Analysis Time	15 samples: <25 min, 96 samples: <140 min
Samples per consumable	15
Sample volume required	1 μ L
Kit stability	4 months
Kit size	Max. 105 samples

¹ Signal-to-noise >3 (single peak)

² Determined using the Genomic DNA ladder as sample

³ Average result from various genomic DNA sample types

⁴ DIN - DNA Integrity Number



Agilent Genomic DNA ScreenTape Assay Quick Guide for 4200 TapeStation System

Storage Conditions

- Reagent vials and the ScreenTape device: 2 – 8 °C (36 – 46 °F).
- Store partially used ScreenTape device upright at 2 – 8 °C (36 – 46 °F) for a maximum of 2 weeks.
- Never freeze ScreenTape device. Discard any accidentally frozen ScreenTape device.

Kit Components

Part Number	Name	Color	Amount
5067-5365	Genomic DNA ScreenTape		7 ScreenTape devices
5067-5366	Genomic DNA Reagents		2 vials
	• Genomic DNA Ladder	●	25 µL
	• Genomic DNA Sample Buffer	●	1350 µL

Limited Use Label License

Some products within this system contain SYBR® Green I, which is licensed from Molecular Probes Inc. for use in research and development only. SYBR® is a registered trademark of Molecular Probes, Inc.

For Research Use Only

Not for use in Diagnostic Procedures.

Additional Material Required for the 4200 TapeStation Instrument

- Loading tips (5067-5598, 1pk or 5067-5599, 10pk)
- 96-well Plates (5042-8502) and 96-well Plate Foil Seal (5067-5154)
- Optical Tube 8x Strip (401428) and Optical Cap 8x Strip (401425)
- Vortex mixer IKA MS3 with adapter

Additional Equipment Required (Not Supplied)

- Volumetric pipette
- Centrifuges for tube strips and well plates

WARNING

Toxic agents

- Refer to product material safety datasheets for further information.
- When working with the ScreenTape assay follow the appropriate safety procedures such as wearing goggles, safety gloves and protective clothing.

CAUTION

Damage to the 4200 TapeStation instrument

- Use only the recommended consumables and reagents with the 4200 TapeStation system.

Essential Measurement Practices


Environmental conditions	<ul style="list-style-type: none">• Optimal operating temperature: 20 °C (68 °F)• Ambient operating temperature: 15 - 30 °C (59 - 86 °F)• Keep reagents during sample preparation at room temperature.
Steps before sample preparation	<ul style="list-style-type: none">• For best results ensure that all kit components are allowed to equilibrate to room temperature for 30 min prior to use.• Vortex mix each vial and briefly spin down.• 'Flick' ScreenTape device to eliminate bubbles in the separation channel, which could interfere with sample loading.
Ladder handling	<ul style="list-style-type: none">• Do not shake or over mix ladder vial, this could result in degradation of the Genomic DNA ladder.
Pipette carefully	<ul style="list-style-type: none">• Always pipette reagents against the side of the well plate or sample tube.• When pipetting small volumes ensure that no liquid remains within the tip.• Ensure that no residual material is left on the outside of the tip.• Care must be taken due to the viscosity of sample buffer.
Mix properly after each pipetting step	<ul style="list-style-type: none">• When adding sample buffer to sample, please ensure that they are mixed correctly.• To achieve this, seal the well plate or cap the sample tubes, centrifuge to collect content at the base, then vortex mix using IKA vortexer and adaptor at 2000 rpm for 1 min.• Improper mixing can lead to quantification errors.
Centrifuge samples before use	<ul style="list-style-type: none">• Briefly centrifuge to ensure all liquid is collected at the bottom of the tubes or well plate and any air bubble is removed.• Apply seal foils to well plates prior to centrifugation and during analysis.• Improper centrifugation of well plates or sample tubes can lead to missing sample data.
Storage after use	<ul style="list-style-type: none">• Reagent vials and the ScreenTape device: 2 – 8 °C (36 – 46 °F).• Store partially used ScreenTape device upright at 2 – 8 °C (36 – 46 °F) for a maximum of 2 weeks.• Never freeze ScreenTape device. Discard any accidentally frozen ScreenTape device.

Ladder considerations

- Ladder is exclusively loaded from location A1 on the tube strip holder.
- The analysis of one ladder per ScreenTape device is required, for 2 – 8 ScreenTape devices a distinct higher ladder volume is prepared.

Agilent Genomic DNA Assay Operating Procedure

- 1 Allow Genomic DNA Reagents (5067-5366) to equilibrate at room temperature for 30 minutes.
- 2 Launch the Agilent 4200 TapeStation Controller Software.
- 3 Flick the Genomic DNA ScreenTape device (5067-5365) and load it into the 4200 TapeStation instrument.
- 4 Place loading tips (5067-5598) into the Agilent 4200 TapeStation instrument.
- 5 Vortex reagents and spin down before use. Take care to not over mix ladder vial.
- 6 Prepare ladder:
 - For 1 – 15 samples: pipette 10 μ L Genomic DNA Sample Buffer (●) and 1 μ L Genomic DNA Ladder (●) at position A1 in a tube strip (401428).
 - For 16 or more samples: pipette 20 μ L Genomic DNA Sample Buffer (●) and 2 μ L Genomic DNA Ladder (●) at position A1 in a tube strip.
- 7 For each sample, pipette 10 μ L Genomic DNA Sample Buffer (●) and 1 μ L DNA sample in a well plate (5042-8502) or a tube strip (401428).
- 8 Apply foil seal (5067-5154) to sample well plate and caps (401425) to tube strips with ladder or sample.
- 9 Mix liquids in sample and ladder vials using the IKA vortex at 2000 rpm for 1 min.
- 10 Spin down to position the sample and ladder at the bottom of the well plate and tube strip.

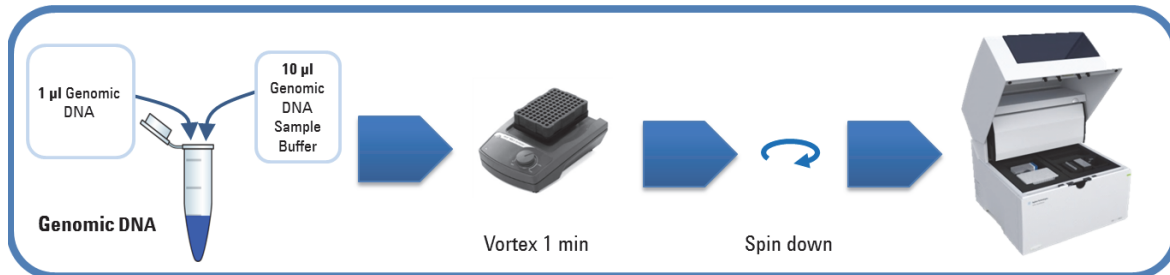

Sample Analysis

- 1 Load samples into the Agilent 4200 TapeStation instrument. Carefully remove caps of tube strips.
- 2 Place ladder in position A1 on tube strip holder in the 4200 TapeStation instrument.
- 3 Select required sample positions on the 4200 TapeStation Controller Software.
- 4 Click **Start**.
- 5 The Agilent TapeStation Analysis Software opens after the run and displays results.

Technical Support and Further Information

For technical support, please visit www.agilent.com/genomics/contact. Visit Agilent Technologies` web site. It offers useful information, support and current developments about the products and technology: www.agilent.com/genomics/tapestation.



G2991-90040

Part Number: G2991-90040

Edition 08/2015

Printed in Germany

© Agilent Technologies, Inc. 2015

Agilent Technologies

Hewlett-Packard-Straße 8

76337 Waldbronn, Germany