



# **SureSelect<sup>QXT</sup> Automated Whole Genome Library Prep for Illumina Multiplexed Sequencing**

**Featuring Transposase-Based Library  
Prep Technology**

**Automated using Agilent NGS  
Workstation Option B**

## **Protocol**

**Version C0, November 2015**

**SureSelect platform manufactured with Agilent  
SurePrint Technology**

**For Research Use Only. Not for use in diagnostic  
procedures.**



**Agilent Technologies**

## Notices

© Agilent Technologies, Inc. 2014, 2015

No part of this manual may be reproduced in any form or by any means (including electronic storage and retrieval or translation into a foreign language) without prior agreement and written consent from Agilent Technologies, Inc. as governed by United States and international copyright laws.

### Manual Part Number

G9682-90010

### Edition

Version C0, November 2015

Printed in USA

Agilent Technologies, Inc.  
5301 Stevens Creek Blvd  
Santa Clara, CA 95051 USA

### Acknowledgement

Oligonucleotide sequences © 2006, 2008, and 2011 Illumina, Inc. All rights reserved. Only for use with the Illumina sequencer systems and associated assays.

### Technical Support

For technical product support, contact your local Agilent Support Services representative.

For US and Canada, call (800) 227-9770 (option 3,4,4). For other countries, find your support center telephone numbers at [www.agilent.com/chem/contactus](http://www.agilent.com/chem/contactus).

Or send an e-mail to:  
[SureSelect.Support@agilent.com](mailto:SureSelect.Support@agilent.com)

## Warranty

**The material contained in this document is provided “as is,” and is subject to being changed, without notice, in future editions. Further, to the maximum extent permitted by applicable law, Agilent disclaims all warranties, either express or implied, with regard to this manual and any information contained herein, including but not limited to the implied warranties of merchantability and fitness for a particular purpose. Agilent shall not be liable for errors or for incidental or consequential damages in connection with the furnishing, use, or performance of this document or of any information contained herein. Should Agilent and the user have a separate written agreement with warranty terms covering the material in this document that conflict with these terms, the warranty terms in the separate agreement shall control.**

## Technology Licenses

The hardware and/or software described in this document are furnished under a license and may be used or copied only in accordance with the terms of such license.

## Restricted Rights Legend

U.S. Government Restricted Rights. Software and technical data rights granted to the federal government include only those rights customarily provided to end user customers. Agilent provides this customary commercial license in Software and technical data pursuant to FAR 12.211 (Technical Data) and 12.212 (Computer Software) and, for the Department of Defense, DFARS 252.227-7015 (Technical Data - Commercial Items) and DFARS 227.7202-3 (Rights in Commercial Computer Software or Computer Software Documentation).

## Notice to Purchaser

This product is provided under an agreement between Bio-Rad Laboratories and Agilent Technologies, Inc., and the manufacture, use, sale or import of this product is subject to US. Pat. No. 6,627,424 and EP Pat. No. 1 283 875 B1, owned by Bio-Rad Laboratories, Inc. Purchase of this product conveys to the buyer the non-transferable right to use the purchased amount of the product and components of the product in PCR (but not real-time PCR) in the Research Field including all Applied Research Fields (including but not limited to forensics, animal testing, and food testing).

## Safety Notices

### CAUTION

A **CAUTION** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in damage to the product or loss of important data. Do not proceed beyond a **CAUTION** notice until the indicated conditions are fully understood and met.

---

### WARNING

A **WARNING** notice denotes a hazard. It calls attention to an operating procedure, practice, or the like that, if not correctly performed or adhered to, could result in personal injury or death. Do not proceed beyond a **WARNING** notice until the indicated conditions are fully understood and met.

---

## In this Guide...

This guide describes an automated protocol for whole genome Illumina paired-end multiplexed library preparation using the SureSelect<sup>QXT</sup> Library Prep system. Sample processing steps are automated using Agilent's NGS Workstation.

### 1 Before You Begin

This chapter contains information that you should read and understand before you start an experiment.

### 2 Using the Agilent NGS Workstation for SureSelect<sup>QXT</sup> Whole Genome Library Prep

This chapter contains an orientation to the Agilent NGS Workstation, an overview of the SureSelect whole genome protocol, and considerations for designing SureSelect experiments for automated processing using the Agilent NGS Workstation.

### 3 Sample Preparation

This chapter describes the steps to prepare gDNA whole genome sequencing libraries.

### 4 Reference

This chapter contains reference information, including component kit contents and index sequences.

## What's New in Version C0

- Updates to custom sequencing primer mixtures for the NextSeq 500 v2 platform (see [Table 30](#) and [Table 31](#) on page 64)
- Update to P5 Index details for NextSeq 500 platform runs using BaseSpace (see [Table 38](#) on page 73)
- Update to Qubit dsDNA Assay Kit nomenclature ([Table 1](#) on page 12 and [step 4](#) on [page 34](#))
- Updates to SureCycler 8800 PCR plate compatibility considerations (see *Caution* on [page 31](#))
- Correction to ordering information for Axygen 96 Deep Well plates (see [Table 2](#) on page 13)
- Revised ordering information for nucleic acid surface decontamination wipes ([Table 2](#) on page 13)

## What's New in Version B1

- Support for sequencing using NextSeq 500 v2 (see [page 61](#) to [page 63](#)).
- Updates to sequencing run setup instructions for 8 bp dual indexes (see [Table 32](#) on page 67).
- Update to Qubit dsDNA Assay Kit ordering information ([Table 1](#) on page 12).

## What's New in Version B0

- Support for new SureSelect<sup>QXT</sup> NSQ Reagent Kits for preparation of DNA libraries for sequencing using Illumina's NextSeq 500 platform. For kit configuration information, see [Table 1](#) on page 12 and [Table 36](#) on page 71. For information on downstream sequencing steps for the NextSeq platform, see "[Step 7. Prepare sequencing samples](#)" on page 61 and see P5 index sequences in [Table 39](#) on page 73.
- Updates to custom sequencing primer preparation instructions for the MiSeq platform ([Table 29](#) on page 63)

- Updates to indexing PCR master mix source plate preparation instructions (changes to order of addition of components, see [page 45](#) to [page 47](#))

# Content

<b>1</b>	<b>Before You Begin</b>	<b>9</b>
	Procedural Notes	10
	Safety Notes	11
	Required Reagents	12
	Required Equipment	13
<b>2</b>	<b>Using the Agilent NGS Workstation for SureSelect<sup>OXT</sup> Whole Genome Library Prep</b>	<b>15</b>
	About the Agilent NGS Workstation	16
	About the Bravo Platform	16
	VWorks Automation Control Software	21
	Overview of the SureSelect <sup>OXT</sup> Whole Genome Library Prep Procedure	27
	Experimental Setup Considerations for Automated Runs	29
	Considerations for Placement of gDNA Samples in 96-well Plates for Automated Processing	30
	Considerations for Equipment Setup	30
	PCR Plate Type Considerations	31
<b>3</b>	<b>Sample Preparation</b>	<b>33</b>
	Step 1. Prepare the genomic DNA samples and Library Prep reagents	34
	Step 2. Fragment and adaptor-tag the genomic DNA samples	36
	Step 3. Amplify adaptor-ligated libraries	43
	Step 4. Purify amplified DNA using AMPure XP beads	53
	Step 5. Assess library DNA quantity and quality using the 2100 Bioanalyzer and High Sensitivity DNA Assay	57
	Step 6. Pool samples for multiplexed sequencing	59
	Step 7. Prepare sequencing samples	61
	Step 8. Set up the sequencing run and trim adaptors from the reads	65

## Contents

<b>4</b>	<b>Reference</b>	<b>69</b>
	Kit Contents	70
	Nucleotide Sequences of SureSelect <sup>QXT</sup> Dual Indexes	72
	Guidelines for Multiplexing with Dual-Indexed Samples	74





# 1 Before You Begin

Procedural Notes	10
Safety Notes	11
Required Reagents	12
Required Equipment	13

Make sure you read and understand the information in this chapter and have the necessary equipment and reagents listed before you start an experiment.

## NOTE

Agilent cannot guarantee the SureSelect kits and cannot provide technical support for the use of non-Agilent protocols or instruments to process samples.



## Procedural Notes

- The SureSelect<sup>QXT</sup> system requires high-quality DNA samples for optimal performance. Use best practices for verifying DNA sample quality before initiating the workflow. For best practice, store diluted DNA solutions at 4°C to avoid repeated freeze-thaw cycles, which may compromise DNA quality.
- Performance of the SureSelect<sup>QXT</sup> library preparation protocol is very sensitive to variations in amounts of DNA sample and other reaction components. It is important to quantify and dilute DNA samples as described on [page 34](#). Carefully measure volumes for all reaction components, and combine components as described in this instruction manual. Use best-practices for liquid handling, including regular pipette calibration, to ensure precise volume measurement.
- Use care in handling the SureSelect QXT Enzyme Mix. After removing the vial from storage at -20°C, keep on ice or in a cold block while in use. Return the vial to storage at -20°C promptly after use.
- Use best-practices to prevent PCR product contamination of samples throughout the workflow:
  - 1 Assign separate pre-PCR and post-PCR pipettors, supplies, and reagents. In particular, never use materials designated to post-PCR segments for the pre-PCR segments of the workflow. For the pre-PCR workflow steps, always use dedicated pre-PCR pipettors with nuclease-free aerosol-resistant tips to pipette dedicated pre-PCR solutions.
  - 2 Maintain clean work areas. Clean pre-PCR surfaces that pose the highest risk of contamination daily using a 10% bleach solution.
  - 3 Wear powder-free gloves. Use good laboratory hygiene, including changing gloves after contact with any potentially-contaminated surfaces.
- Possible stopping points, where samples may be stored at -20°C, are marked in the protocol. Do not subject the samples to multiple freeze/thaw cycles.
- To prevent contamination of reagents by nucleases, always wear powder-free laboratory gloves and use dedicated solutions and pipettors with nuclease-free aerosol-resistant tips.
- In general, follow Biosafety Level 1 (BL1) safety rules.

## Safety Notes

### CAUTION

- Wear appropriate personal protective equipment (PPE) when working in the laboratory.
-

## Required Reagents

**Table 1** Required Reagents for SureSelect<sup>OXT</sup> Whole Genome Library Prep

Description	Vendor and part number
SureSelect <sup>OXT</sup> Library Prep Kit	Agilent
Illumina HiSeq or MiSeq platform (ILM), 96 Samples	p/n G9682B
Illumina NextSeq platform (NSQ), 96 Samples	p/n G9684B
Agencourt AMPure XP Kit	Beckman Coulter Genomics
5 mL	p/n A63880
60 mL	p/n A63881
450 mL	p/n A63882
1X Low TE Buffer (10 mM Tris-HCl, pH 8.0, 0.1 mM EDTA)	Life Technologies p/n 12090-015, or equivalent
100% Ethanol, molecular biology grade	Sigma-Aldrich p/n E7023
Qubit dsDNA HS Assay Kit <i>or</i>	Life Technologies p/n Q32851
Qubit dsDNA BR Assay Kit	Life Technologies
100 assays	p/n Q32850
500 assays	p/n Q32853
Nuclease-free Water (not DEPC-treated)	Ambion Cat #AM9930

## Required Equipment

**Table 2** Required Equipment for SureSelect<sup>QXT</sup> Whole Genome Library Prep

Description	Vendor and part number
Agilent NGS Workstation Option B, with VWorks software version 11.3.0.1195	Agilent p/n G5522A Contact Agilent Automation Solutions for more information: Customerservice.automation@agilent.com
Robotic Pipetting Tips (Sterile, Filtered, 250 µL)	Agilent p/n 19477-022
Clear Peelable Seal plate seals (for use with the PlateLoc Thermal Plate Sealer)	Agilent p/n 16985-001
Thermal cycler and accessories	SureCycler 8800 Thermal Cycler (Agilent p/n G8810A), 96 well plate module (Agilent p/n G8810A) and compression mats (Agilent p/n 410187) or equivalent
PCR plates compatible with selected Thermal Cycler, e.g. Agilent semi-skirted PCR plate for the SureCycler 8800 Thermal Cycler When selecting plates for another thermal cycler, see <a href="#">Table 6</a> on page 31 for the list of PCR plates supported in automation protocols	Agilent p/n 401334
Eppendorf twin.tec full-skirted 96-well PCR plates	Eppendorf p/n 951020401 or 951020619
Thermo Scientific Reservoirs	Thermo Scientific p/n 1064156
Nunc DeepWell Plates, sterile, 1.3-mL well volume	Thermo Scientific p/n 260251
Axygen 96 Deep Well Plate, 2 mL, Square Well (waste reservoirs; working volume 2.2 mL)	Axygen p/n P-2ML-SQ-C E & K Scientific p/n EK-2440
Nucleic acid surface decontamination wipes	DNA Away Surface Decontaminant Wipes, Thermo Scientific p/n 7008, or equivalent
<b>DNA Analysis Platform and Consumables</b>	
Agilent 2100 Bioanalyzer Laptop Bundle	Agilent p/n G2943CA
Agilent 2100 Bioanalyzer Electrophoresis Set	Agilent p/n G2947CA
Agilent DNA 1000 Kit	Agilent p/n 5067-1504
Agilent High Sensitivity DNA Kit	Agilent p/n 5067-4626

## 1 Before You Begin

### Required Equipment

**Table 2** Required Equipment for SureSelect<sup>OXT</sup> Whole Genome Library Prep

Description	Vendor and part number
Qubit Fluorometer	Life Technologies p/n Q32857
Qubit Assay Tubes	Life Technologies p/n Q32856
DNA LoBind Tubes, 1.5-mL PCR clean, 250 pieces	Eppendorf p/n 022431021 or equivalent
Centrifuge	Eppendorf Centrifuge model 5804 or equivalent
Plate or strip tube centrifuge	Labnet International MPS1000 Mini Plate Spinner p/n C1000 (requires adapter, p/n C1000-ADAPT, for use with strip tubes) or equivalent
Multichannel pipette	Pipetman or equivalent
P10, P20, P200 and P1000 pipettes	Pipetman P10, P20, P200, P1000 or equivalent
Vortex mixer	
Ice bucket	
Powder-free gloves	
Sterile, nuclease-free aerosol barrier pipette tips	



## 2 Using the Agilent NGS Workstation for SureSelect<sup>QXT</sup> Whole Genome Library Prep

About the Agilent NGS Workstation 16

Overview of the SureSelect<sup>QXT</sup> Whole Genome Library Prep  
Procedure 27

Experimental Setup Considerations for Automated Runs 29

This chapter contains an orientation to the Agilent NGS Workstation, an overview of the SureSelect<sup>QXT</sup> whole genome library preparation protocol, and considerations for designing SureSelect<sup>QXT</sup> experiments for automated processing using the Agilent NGS Workstation.



## About the Agilent NGS Workstation

### About the Bravo Platform


The Bravo platform is a versatile liquid handler with a nine plate-location platform deck, suitable for handling 96-well, 384-well, and 1536-well plates. The Bravo platform is controlled by the VWorks Automation Control software. Fitted with a choice of seven interchangeable fixed-tip or disposable-tip pipette heads, it accurately dispenses fluids from 0.1  $\mu\text{L}$  to 250  $\mu\text{L}$ .

#### CAUTION

Before you begin, make sure that you have read and understand operating, maintenance and safety instructions for using your Bravo platform. Refer to the *Bravo Platform User Guide* (G5409-90006) and the *VWorks Software User Guide* (G5415-90063).

#### Bravo Platform Deck

The protocols in the following sections include instructions for placing plates and reagent reservoirs on specific Bravo deck locations. Use [Figure 1](#) to familiarize yourself with the location numbering convention on the Bravo platform deck.

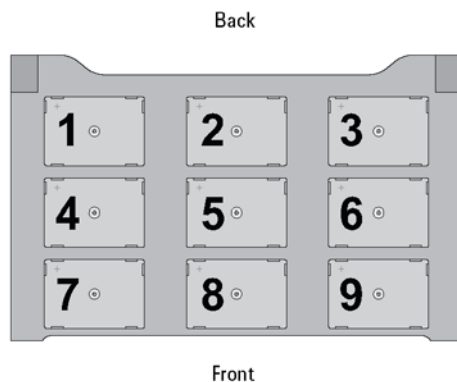

**Figure 1** Bravo platform deck



### Setting the Temperature of Bravo Deck Heat Blocks


Bravo deck positions 4 and 6 are equipped with Inheco heat blocks, used to incubate sample plates at defined temperatures during the run. Runs that include high- (85°C) or low- (4°C) temperature incubation steps may be expedited by pre-setting the temperature of the affected block before starting the run.

Bravo deck heat block temperatures may be changed using the Inheco Multi TEC Control device touchscreen as described in the steps below. See [Table 3](#) for designations of the heat block-containing Bravo deck positions on the Multi TEC control device.

**Table 3** Inheco Multi TEC Control touchscreen designations


Bravo Deck Position	Designation on Inheco Multi TEC Control Screen
4	CPAC 2 1
6	CPAC 2 2

- Using the arrow buttons, select the appropriate block (CPAC 2 block 1 or CPAC 2 block 2).

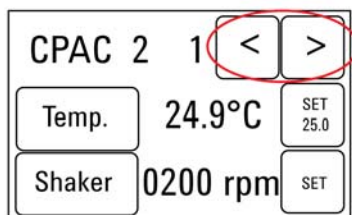


## 2 Using the Agilent NGS Workstation for SureSelect<sup>QXT</sup> Whole Genome Library Prep About the Bravo Platform


- 2 To set the temperature of the selected block, press the SET button.

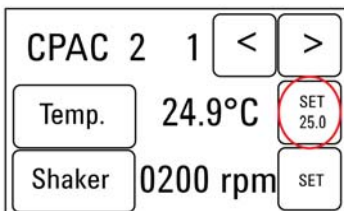

- 3 Using the numeral pad, enter the desired temperature. The entered temperature appears in the top, left rectangle. Once the correct temperature is displayed, press the rectangle to enter the temperature.

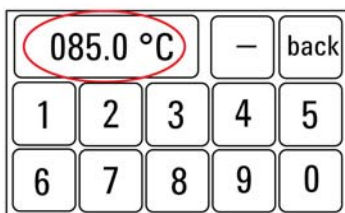

- 4 Press the Temp button until the new temperature is displayed on the SET button and until the Temp button is darkened, indicating that the selected heat block is heating or cooling to the new temperature setting. The current temperature of the block is indicated in the center of the display.

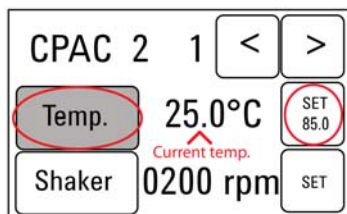

### Setting the Temperature of Bravo Deck Position 9 Using the ThermoCube Device

Bravo deck position 9 is equipped with a ThermoCube thermoelectric temperature control system, used to incubate components at a defined temperature during the run. During protocols that require temperature control at position 9, you will be instructed to start and set the temperature of the ThermoCube device before starting the run.

ThermoCube temperature settings are modified using the control panel (LCD display screen and four input buttons) on the front panel of the device using the following steps.

- 1 Turn on the ThermoCube and wait for the LCD screen to display **TEMP.**
- 2 Press the **UP** or **DOWN** button to change **SET TEMP 1** to the required set point.
- 3 Press the **START** button.


The ThermoCube then initiates temperature control of Bravo deck position 9 at the displayed set point.

## 2 Using the Agilent NGS Workstation for SureSelect<sup>0XT</sup> Whole Genome Library Prep About the Bravo Platform

### Using the Labware MiniHub

The protocols in the following sections include instructions for placing plates or reservoirs at specific Labware MiniHub positions. Use [Figure 2](#) to familiarize yourself with the required orientations loading plates in the Labware MiniHub for use in SureSelect automation protocols.

For Thermo Scientific reservoirs, place the notched corner facing the center of the hub.

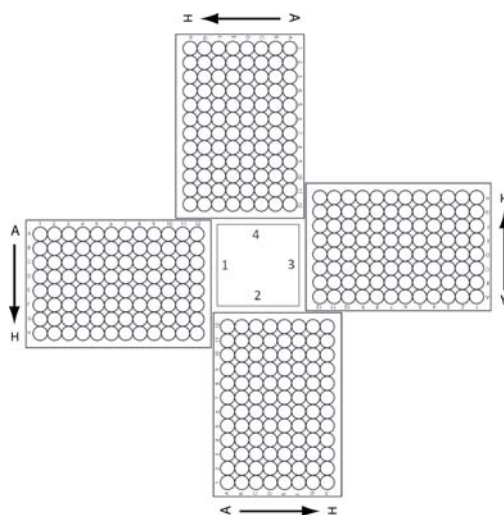

**Figure 2** Agilent Labware MiniHub plate orientation.

## VWorks Automation Control Software

VWorks software, included with your Agilent NGS Workstation, allows you to control the robot and integrated devices using a PC. The Agilent NGS Workstation is preloaded with VWorks software containing all of the necessary SureSelect system liquid handling protocols. General instructions for starting up the VWorks software and the included protocols is provided below. Each time a specific VWorks protocol is used in the SureSelect procedure, any settings required for that protocol are included in the relevant section of this manual.

### NOTE

The instructions in this manual are compatible with VWorks software version 11.3.0.1195, including SureSelect<sup>QXT</sup> automation protocols version 1.0.

If you have questions about VWorks version compatibility, please contact [service.automation@agilent.com](mailto:service.automation@agilent.com).

### Logging in to the VWorks software

- 1 Double-click the VWorks icon or the SureSelectQXT\_ILM\_v1.0.VWForm shortcut on the Windows desktop to start the VWorks software.
- 2 If User Authentication dialog is not visible, click **Log in** on the VWorks window toolbar.
- 3 In the User Authentication dialog, type your VWorks user name and password, and click **OK**. (If no user account is set up, contact the administrator.)


### VWorks protocol and runset files

VWorks software uses two file types for automation runs, .pro (protocol) files and .rst (runset) files. Runset files are used for automated procedures in which the workstation uses more than one automation protocol during the run.

## 2 Using the Agilent NGS Workstation for SureSelect<sup>QXT</sup> Whole Genome Library Prep VWorks Automation Control Software

### Using the SureSelect<sup>QXT</sup>\_ILM\_v1.0.VWForm to setup and start a run

Use the VWorks form SureSelect<sup>QXT</sup>\_ILM\_v1.0.VWForm, shown below, to set up and start each SureSelect automation protocol or runset.



**SureSelect<sup>QXT</sup>**  
Transposase Library Prep  
for Illumina sequencers

**Parameters**

1) Select Protocol to Run

AMPureXP Case:

2) Select PCR Plate/labware for Thermal Cycling

3) Select Number of Columns of Samples

4) Click button below to Display Initial Workstation Setup

5) Load labware according to Workstation Setup -->

**Controls**

Once you have loaded labware according to Workstation Setup on right, click "Run Selected Protocol" to start run.

Elapsed Time: 00:00:00

**Information**

Currently Running Protocol:

**Advanced Settings**

TESTING ONLY: Reduces all incubation times

**Workstation Setup**

MiniHub	MiniHub Cassette 1	MiniHub Cassette 2	MiniHub Cassette 3	MiniHub Cassette 4
Shelf 5				
Shelf 4				
Shelf 3				
Shelf 2				
Shelf 1				


**Bravo Deck**

<Position 1>	<Position 2>	<Position 3>
<Pos 4: Peltier>	<Pos 5: Shaker>	<Pos 6: Peltier>
<Pos 7: Magnetic>	<Position 8>	<Pos 9: Chiller>

**BenchCel**

BenchCel Stacker 1	BenchCel Stacker 2	BenchCel Stacker 3	BenchCel Stacker 4

- 1 Open the form using the SureSelect<sup>QXT</sup>\_ILM\_v1.0.VWForm shortcut on your desktop.
- 2 Use the form drop-down menus to select the appropriate SureSelect workflow step and number of columns of samples for the run.
- 3 Once all run parameters have been specified on the form, click **Display Initial Workstation Setup**.

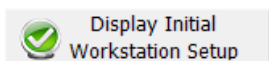

- 4 The Workstation Setup region of the form will then display the required placement of reaction components and labware on the NGS Workstation for the specified run parameters.



**SureSelect<sup>QXT</sup>**  
Transposase Library Prep  
for Illumina sequencers

**Parameters**

1) Select Protocol to Run

LibraryPrep\_QXT\_ILM\_v1.0.rst

AMPureXP Case: Not Applicable

2) Select PCR Plate labware for Thermal Cycling

96 Agilent Semi-skirted PCR in Red Alum Insert

3) Select Number of Columns of Samples

12

4) Click button below to Display Initial Workstation Setup

Display Initial Workstation Setup

Clear Workstation Setup Display

5) Load labware according to Workstation Setup -->

**Controls**

Once you have loaded labware according to Workstation Setup on right, click "Run Selected Protocol" to start run.

Run Selected Protocol

Pause

Initialize all devices

Full Screen

Gantt Chart

Elapsed Time: 00:00:00

Reset All Form Selections to Defaults

**Information**

Currently Running Protocol:

**Advanced Settings**

TESTING ONLY: Reduces all incubation times

**Workstation Setup**

MiniHub

	MiniHub Cassette 1	MiniHub Cassette 2	MiniHub Cassette 3	MiniHub Cassette 4
Shelf 5		Empty Nunc DeepWell Plate		
Shelf 4				Stop Solution (twin.tec)
Shelf 3		Empty Eppendorf twin.tec Plate		
Shelf 2	New Tip Box	Nuclease-free Water Reservoir	AmpureXP Beads in Nunc DeepWell	
Shelf 1	Empty Tip Box	70% Ethanol Reservoir		Empty Tip Box


Bravo Deck

<Position 1> Waste Reservoir (Axygen 96DW)	<Position 2>	<Position 3>
<Pos 4: Peltier> 52°C Red Insert	<Pos 5: Shaker>	<Pos 6: Peltier> 4°C Nunc MasterMix Plate (Col 1)
<Pos 7: Magnetic> DNA Plate (twin.tec)	<Position 8>	<Pos 9: Chiller> 0°C Empty Eppendorf twin.tec Plate on Red Insert

BenchCel

BenchCel Stacker 1	BenchCel Stacker 2	BenchCel Stacker 3	BenchCel Stacker 4
8 Tip Boxes	Empty	Empty	Empty


- 5 After verifying that the NGS Workstation has been set up correctly, click **Run Selected Protocol**.

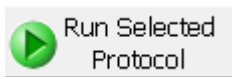

### Error messages encountered at start of run


After starting the run, you may see the error messages displayed below. When encountered, make the indicated selections and proceed with the run. Encountering either or both of these error messages is not indicative of a problem with the NGS workstation or your run setup.

- 1 If you encounter the G-axis error message shown below, select **Ignore and Continue, leaving device in current state**.

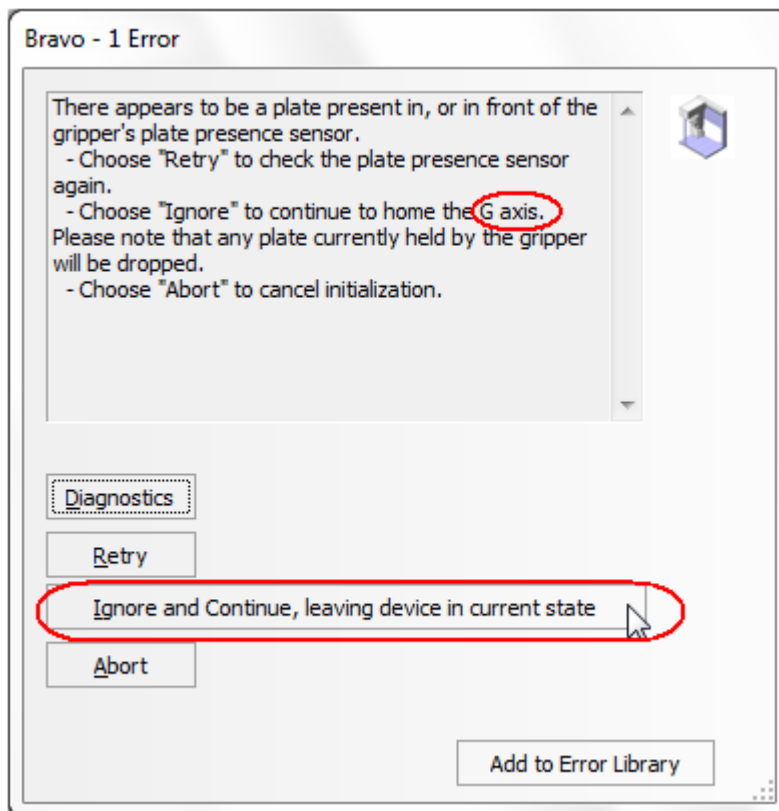



- 2 If you encounter the W-axis error message shown below, select **Retry**.

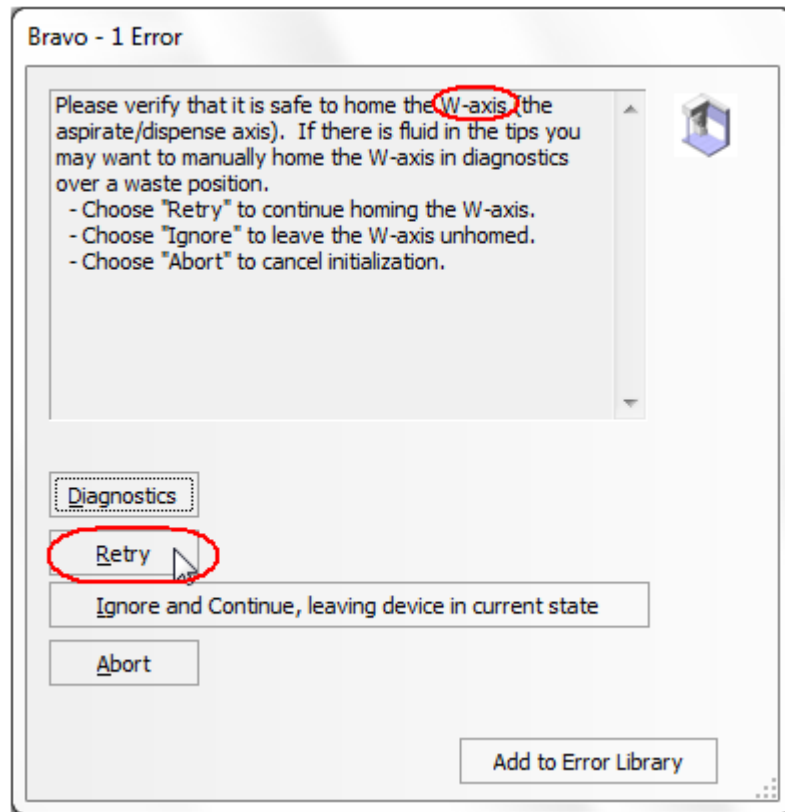

### Verifying the Simulation setting

VWorks software may be run in simulation mode, during which commands entered on screen are not completed by the NGS workstation. If workstation devices do not respond when you start a run, verify the simulation mode status in VWorks using the following steps.

- 1 Verify that **Simulation is off** is displayed on the status indicator (accessible by clicking **View > Control Toolbar**).




- 2 If the indicator displays **Simulation is on**, click the status indicator button to turn off the simulation mode.

#### NOTE

If you cannot see the toolbar above the SureSelect\_XT\_Illumina VWorks form, click the **Full Screen** button to exit full screen mode. If the toolbar is still not visible, right-click on the form and then select **Control Toolbar** from the menu.

### Finishing a protocol or runset

The window below appears when each run is complete. Click **Yes** to release the BenchCel racks to allow removal of components used in the current run in preparation for the next .pro or .rst run.

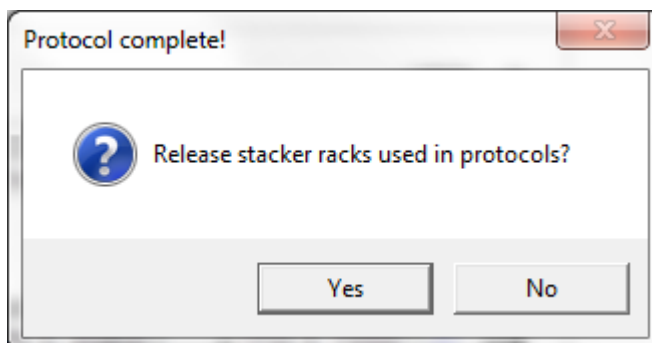

## Overview of the SureSelect<sup>QXT</sup> Whole Genome Library Prep Procedure


Figure 3 summarizes the SureSelect<sup>QXT</sup> whole genome library prep workflow for samples to be sequenced using the Illumina paired-read sequencing platform. Each sample to be sequenced requires an individual library preparation reaction. The samples are then tagged by PCR with dual index sequences. Depending on the experimental design, up to 96 samples can be pooled and sequenced in a single lane using the dual index tags that are provided with SureSelect<sup>QXT</sup> Library Prep kits.

Table 4 summarizes how the VWorks protocols are integrated into the SureSelect<sup>QXT</sup> workflow. See the [Sample Preparation](#) chapter for complete instructions for use of the VWorks protocols for sample processing.

## 2 Using the Agilent NGS Workstation for SureSelect<sup>QXT</sup> Whole Genome Library Prep

Overview of the SureSelect<sup>QXT</sup> Whole Genome Library Prep Procedure

### SureSelect<sup>QXT</sup> Whole Genome NGS Workflow

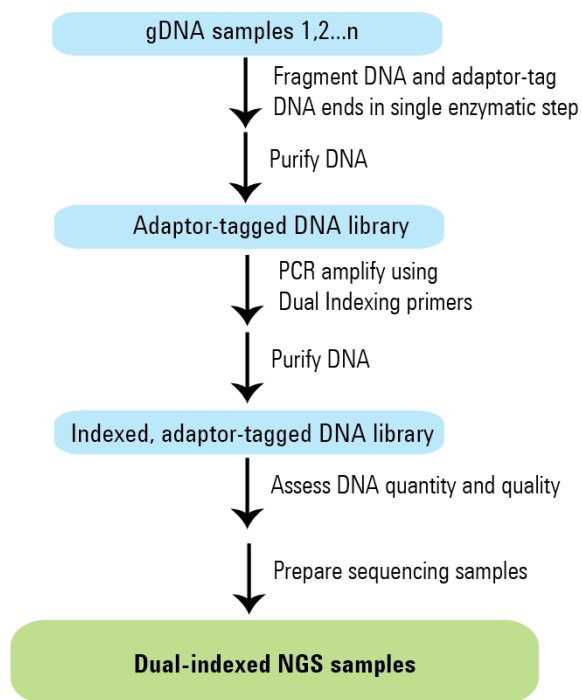

**Figure 3** Overall sequencing sample preparation workflow.

**Table 4** Overview of VWorks protocols and runsets used for SureSelect<sup>QXT</sup> Whole Genome Library Prep

Workflow Step	VWorks Protocol Used for Agilent NGS Workstation automation
Prepare fragmented and adaptor-tagged DNA	LibraryPrep_QXT_ILM_v1.0.rst
Purify adaptor-tagged DNA using AMPure XP beads	
Amplify adaptor-tagged DNA and add index tags	WholeGenomePCR_QXT_ILM_v1.0.pro
Purify indexed DNA using AMPure XP beads	AMPureXP_QXT_ILM_v1.0.pro:Whole Genome PCR

## Experimental Setup Considerations for Automated Runs

Agilent SureSelect Automated Library Prep runs may include 1, 2, 3, 4, 6, or 12 columns (equivalent to 8, 16, 24, 32, 48, or 96 wells) of gDNA samples to be prepared for sequencing on the Illumina platform. Plan your experiments using complete columns of samples.

**Table 5** Columns to Samples Equivalency

Number of Columns Processed	Total Number of Samples Processed
1	8
2	16
3	24
4	32
6	48
12	96

The number of columns or samples that may be processed using the supplied reagents will depend on the experimental design. For greatest efficiency of reagent use, plan experiments using at least 3 columns per run. Each 96-reaction kit contains sufficient reagents for 96 reactions configured as 4 runs of 3 columns of samples per run.

## Considerations for Placement of gDNA Samples in 96-well Plates for Automated Processing

- The Agilent NGS Workstation processes samples column-wise beginning at column 1. gDNA samples should be loaded into 96-well plates column-wise, in well order A1 to H1, then A2 to H2, ending with A12 to H12. When processing partial runs with <12 sample columns, do not leave empty columns between sample columns; always load the plate using the left-most column that is available.
- Dual index assignments for the DNA samples can affect sample placement decisions at the beginning of the workflow. For example, all samples on the same row of the DNA sample plate must be assigned to the same P5 indexing primer during sample indexing (see [Figure 3](#)). It is important to review and understand the guidelines for assignment of dual indexing primers on [page 45](#) while planning sample placement for the run to ensure that the indexing design is compatible with the initial DNA sample placement.


## Considerations for Equipment Setup

- Some workflow steps require the rapid transfer of sample plates between the Bravo deck and a thermal cycler. Locate your thermal cycler in close proximity to the Agilent NGS Workstation to allow rapid and efficient plate transfer.
- Several workflow steps require that the sample plate be sealed using the PlateLoc thermal microplate sealer included with the Agilent NGS Workstation, and then centrifuged to collect any dispersed liquid. To maximize efficiency, locate the centrifuge in close proximity to the Agilent NGS Workstation.

## PCR Plate Type Considerations

Automation protocols include liquid-handling steps in which reagents are dispensed to PCR plates in preparation for transfer to a thermal cycler. For these steps you must specify the PCR plate type to be used on the SureSelectQXT\_ILM\_v1.0.VWForm to allow correct configuration of the liquid handling components for the PCR plate type. Before you begin the automation protocol, make sure that you are using a supported PCR plate type. The PCR plate type to be used in the protocol is specified using the menu below. Vendor and part number information is provided for the supported plate types in [Table 6](#).

### 2) Select PCR Plate labware for Thermal Cycling

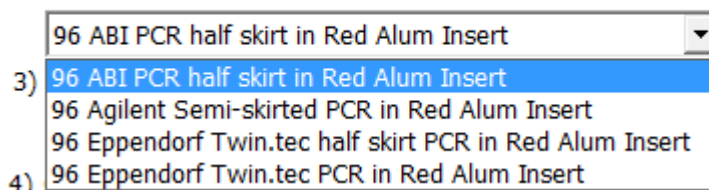

**Table 6** Ordering information for supported PCR plates

Description in VWorks menu	Vendor and part number
96 ABI PCR half-skirted plates (MicroAmp Optical plates)	Life Technologies p/n N8010560
96 Agilent semi-skirted PCR plate	Agilent p/n 401334
96 Eppendorf Twin.tec half-skirted PCR plates	Eppendorf p/n 951020303
96 Eppendorf Twin.tec PCR plates (full-skirted)	Eppendorf p/n 951020401

### CAUTION

The plates listed in [Table 6](#) are compatible with the Agilent NGS Bravo and associated VWorks automation protocols, designed to support use of various thermal cyclers.

Accordingly, some plates listed in [Table 6](#) are not compatible with the recommended SureCycler 8800 Thermal Cycler. When using the SureCycler 8800 Thermal Cycler in the SureSelect automation workflow, use 96 Agilent semi-skirted PCR plates.

When using a different thermal cycler in the workflow, be sure to select a PCR plate that is compatible with your thermal cycler and that is listed in [Table 6](#).

## **2 Using the Agilent NGS Workstation for SureSelect<sup>0XT</sup> Whole Genome Library Prep** **PCR Plate Type Considerations**





## 3 Sample Preparation

- Step 1. Prepare the genomic DNA samples and Library Prep reagents 34
- Step 2. Fragment and adaptor-tag the genomic DNA samples 36
- Step 3. Amplify adaptor-ligated libraries 43
- Step 4. Purify amplified DNA using AMPure XP beads 53
- Step 5. Assess library DNA quantity and quality using the 2100 Bioanalyzer and High Sensitivity DNA Assay 57
- Step 6. Pool samples for multiplexed sequencing 59
- Step 7. Prepare sequencing samples 61
- Step 8. Set up the sequencing run and trim adaptors from the reads 65

This section contains instructions for whole genome library preparation from gDNA specific to the Illumina paired-read sequencing platform and to automated processing using the Agilent NGS Workstation.



### 3 Sample Preparation

#### Step 1. Prepare the genomic DNA samples and Library Prep reagents

## Step 1. Prepare the genomic DNA samples and Library Prep reagents

It is important to have all materials prepared in advance of use in the SureSelect<sup>QXT</sup> automated Library Prep protocol. In this step, the gDNA is carefully quantified and dispensed into the sample plate. Additional reagents that require modification or temperature equilibration before use are also prepared in this step.

- 1 Remove the DMSO vial from the SureSelect QXT Library Prep Kit Box 2 in  $-20^{\circ}\text{C}$  storage. Leave the DMSO vial at room temperature in preparation for use on [page 45](#).
- 2 Prepare reagents for the purification protocols on [page 37](#) and [page 53](#).
  - a Transfer the AMPure XP beads to room temperature. The beads should be held at room temperature for at least 30 minutes before use. *Do not freeze the beads at any time.*
  - b Prepare 100 mL of fresh 70% ethanol for use in the purification steps. The 70% ethanol may be used for multiple steps done on the same day, when stored in a sealed container.
- 3 Obtain the bottle of SureSelect QXT Stop Solution from SureSelect QXT Library Prep Box 1 (stored at room temperature). Verify that the SureSelect QXT Stop Solution contains 25% ethanol, by referring to the container label and the instructions below.

Before the first use of a fresh container, add 1.5 mL of ethanol to the provided bottle containing 4.5 mL of stop solution, for a final ethanol concentration of 25%. Seal the bottle then vortex well to mix. After adding the ethanol, be sure to mark the label for reference by later users.

Keep the prepared 1X SureSelect QXT Stop Solution at room temperature, tightly sealed, until it is used on [page 38](#).
- 4 Quantify and dilute gDNA samples using two serial fluorometric assays:
  - a Use the Qubit dsDNA BR Assay or Qubit dsDNA HS Assay to determine the initial concentration of each gDNA sample. Follow the manufacturer's instructions for the specific assay kit and the Qubit

## Step 1. Prepare the genomic DNA samples and Library Prep reagents

instrument. This step is critical for successful preparation of input DNA at the required concentration to ensure optimal fragmentation.

- b** Dilute each gDNA sample with nuclease-free water to a final concentration of 100 ng/μL in a 1.5-mL LoBind tube.
  - c** Carefully measure the DNA concentration of each of the 100 ng/μL dilutions using a second Qubit dsDNA BR or HS Assay.
  - d** Adjust each gDNA sample with nuclease-free water to a final concentration of 10 ng/μL in a 1.5-mL LoBind tube.
- 5** Transfer 5 μL of the 10 ng/μL-DNA samples into the wells of a 96-well Eppendorf plate, column-wise, for processing on the Agilent NGS Workstation, in well order A1 to H1, then A2 to H2, ending with A12 to H12.

**NOTE**

SureSelect Automated Library Prep runs may include 1, 2, 3, 4, 6, or 12 columns of the plate. See [Experimental Setup Considerations for Automated Runs](#) on [page 29](#) for additional sample placement considerations.

- 6** Seal the plate using the PlateLoc Thermal Microplate Sealer, with sealing settings of 165°C and 1.0 sec.
- 7** Centrifuge the plate for 30 seconds to drive the well contents off the walls and plate seal and to remove air bubbles.

Store the sample plate on ice until it is used on [page 41](#).

### 3 Sample Preparation

#### Step 2. Fragment and adaptor-tag the genomic DNA samples

## Step 2. Fragment and adaptor-tag the genomic DNA samples

In this step, automation runset LibraryPrep\_QXT\_ILM\_v1.0.rst is used to enzymatically fragment the gDNA and to add adaptors to ends of the fragments in a single reaction. After fragmentation and tagging, the Agilent NGS Workstation purifies the prepared DNA using AMPure XP beads.

This step uses the SureSelect<sup>QXT</sup> Reagent Kit components listed in [Table 7](#) in addition to reagents prepared for use on [page 34](#) to [page 35](#).

**Table 7** Reagents for DNA fragmentation and adaptor-tagging

Kit Component	Storage Location	Where Used
SureSelect QXT Buffer	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 36</a>
SureSelect QXT Enzyme Mix ILM	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 36</a>

### Prepare the workstation

- 1 Clear the Labware MiniHub and BenchCel of all plates and tip boxes.
- 2 Gently wipe down the Labware MiniHub, Bravo decks, and BenchCel with a DNA Away decontamination wipe.
- 3 Pre-set the temperature of Bravo deck position 4 to 52°C and position 6 to 4°C using the Inheco Multi TEC control touchscreen, as described in [Setting the Temperature of Bravo Deck Heat Blocks](#). On the control touchscreen, Bravo deck position 4 corresponds to CPAC 2, position 1 and Bravo deck position 6 corresponds to CPAC 2, position 2.
- 4 Turn on the ThermoCube, set to 0°C, at position 9 of the Bravo deck. Be sure that the chiller reservoir contains at least 300 mL of 25% ethanol.
- 5 Place red PCR plate inserts at Bravo deck positions 4 and 9.

## Step 2. Fragment and adaptor-tag the genomic DNA samples

- 6 Load tip boxes for the run in the BenchCel Microplate Handling Workstation according to [Table 8](#).

**Table 8** Initial BenchCel configuration for LibraryPrep\_QXT\_ILM\_v1.0.rst

No. of Columns Processed	Rack 1	Rack 2	Rack 3	Rack 4
1	1 Tip box	Empty	Empty	Empty
2	2 Tip boxes	Empty	Empty	Empty
3	2 Tip boxes	Empty	Empty	Empty
4	3 Tip boxes	Empty	Empty	Empty
6	4 Tip boxes	Empty	Empty	Empty
12	8 Tip boxes	Empty	Empty	Empty

- 7 Load the workstation MiniHub with the empty plates and other labware components for the run, using the positions shown in the Workstation Setup region of the VWorks Form. Use the plate orientations shown in [Figure 2](#) on page 20.

### Prepare the purification reagents

- 8 Verify that the AMPure XP bead suspension is at room temperature. *Do not freeze the beads at any time.*
- 9 Mix the bead suspension well so that the reagent appears homogeneous and consistent in color.
- 10 Prepare a Nunc DeepWell source plate for the beads by adding 55  $\mu$ L of homogeneous AMPure XP beads per well, for each well to be processed. Place the bead source plate on shelf 2 of cassette 3 of the workstation MiniHub.
- 11 Prepare a Thermo Scientific reservoir containing 15 mL of nuclease-free water. Place the water reservoir on shelf 2 of cassette 2 of the workstation MiniHub.
- 12 Prepare a separate Thermo Scientific reservoir containing 45 mL of freshly-prepared 70% ethanol. Place the ethanol reservoir on shelf 1 of cassette 2 of the workstation MiniHub.

### 3 Sample Preparation

#### Step 2. Fragment and adaptor-tag the genomic DNA samples

##### Prepare the Library Prep Master Mix and Stop Solution source plates

**13** Prepare the Stop Solution source plate using an Eppendorf twin.tec full-skirted PCR plate. Add 35  $\mu\text{L}$  of 1X SureSelect QXT Stop Solution per well, for each well to be processed. Place the source plate on shelf 4 of cassette 4 of the workstation MiniHub.

**14** Before use, vortex the SureSelect QXT Buffer and SureSelect QXT Enzyme Mix ILM tubes vigorously at high speed.

These components are in liquid form when removed from  $-20^{\circ}\text{C}$  storage and should be returned to  $-20^{\circ}\text{C}$  storage promptly after use.

#### CAUTION

Minor variations in volumes of the solutions combined in [step 15](#) below may result in DNA fragment size variation.

The SureSelect QXT Buffer and Enzyme Mix solutions are highly viscous. Thorough mixing of the reagents is critical for optimal performance.

**15** Prepare the appropriate volume of Library Prep Master Mix, according to [Table 9](#). Mix well by vortexing for 20 seconds and then keep on ice.

**Table 9** Preparation of Library Prep Master Mix

SureSelect <sup>QXT</sup> Reagent	Volume for 1 Library	Volume for 1 Column	Volume for 2 Columns	Volume for 3 Columns	Volume for 4 Columns	Volume for 6 Columns	Volume for 12 Columns
SureSelect QXT Buffer	17.0 $\mu\text{L}$	216.8 $\mu\text{L}$	361.3 $\mu\text{L}$	505.8 $\mu\text{L}$	650.3 $\mu\text{L}$	939.3 $\mu\text{L}$	1878.5 $\mu\text{L}$
Nuclease-free water	1.0 $\mu\text{L}$	12.8 $\mu\text{L}$	21.3 $\mu\text{L}$	29.8 $\mu\text{L}$	38.3 $\mu\text{L}$	55.3 $\mu\text{L}$	110.5 $\mu\text{L}$
SureSelect QXT Enzyme Mix ILM	1.0 $\mu\text{L}$	12.8 $\mu\text{L}$	21.3 $\mu\text{L}$	29.8 $\mu\text{L}$	38.3 $\mu\text{L}$	55.3 $\mu\text{L}$	110.5 $\mu\text{L}$
<b>Total Volume</b>	<b>19 <math>\mu\text{L}</math></b>	<b>242.3 <math>\mu\text{L}</math></b>	<b>403.8 <math>\mu\text{L}</math></b>	<b>565.3 <math>\mu\text{L}</math></b>	<b>726.8 <math>\mu\text{L}</math></b>	<b>1049.8 <math>\mu\text{L}</math></b>	<b>2099.5 <math>\mu\text{L}</math></b>


**16** Prepare the Library Prep master mix source plate using a Nunc DeepWell plate, containing the mixture from [step 15](#). Add the volume indicated in [Table 10](#) to all wells of column 1 of the Nunc DeepWell

Step 2. Fragment and adaptor-tag the genomic DNA samples

plate. Keep the master mix on ice during the aliquoting steps. The final configuration of the master mix source plate is shown in Figure 4.

**Table 10** Preparation of the Master Mix Source Plate for LibraryPrep\_QXT\_ILM\_v1.0.rst

Master Mix Solution	Position on Source Plate	Volume of Master Mix added per Well of Nunc Deep Well Source Plate					
		1-Column Runs	2-Column Runs	3-Column Runs	4-Column Runs	6-Column Runs	12-Column Runs
Library Prep Master Mix	Column 1 (A1-H1)	27.9 µL	48.1 µL	68.3 µL	88.5 µL	128.8 µL	260.1 µL

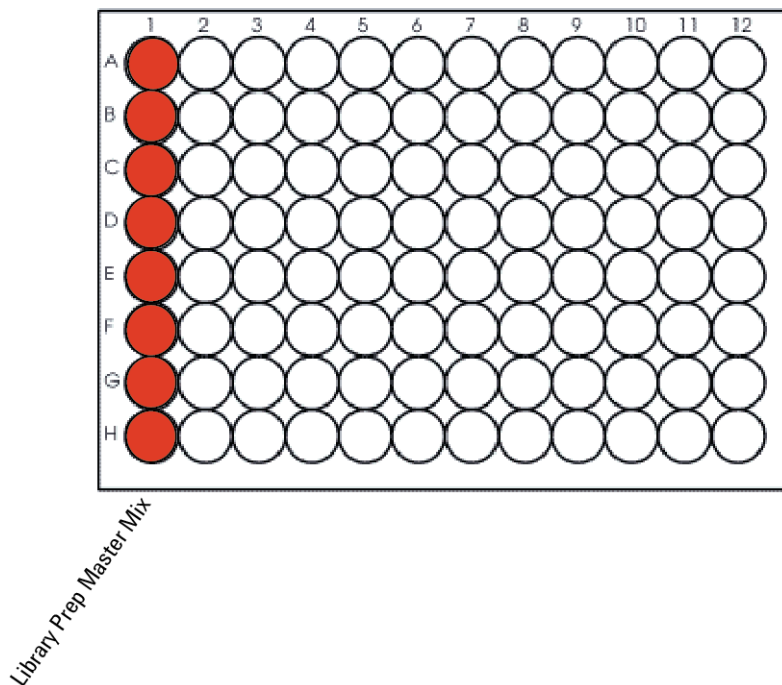

**Figure 4** Configuration of the master mix source plate for LibraryPrep\_QXT\_ILM\_v1.0.rst

### 3 Sample Preparation

#### Step 2. Fragment and adaptor-tag the genomic DNA samples

**17** Seal the master mix source plate using the PlateLoc Thermal Microplate Sealer, with sealing settings of 165°C and 1.0 sec.

**18** Centrifuge the plate for 30 seconds to drive the well contents off the walls and plate seal and to eliminate any bubbles. Keep the master mix source plate on ice.

#### NOTE

The presence of bubbles in source plate solutions may cause inaccurate volume transfer by the Bravo liquid handling platform. Ensure that the source plate is sealed and centrifuged prior to use in a run.

#### Load the Agilent NGS Workstation

**19** Verify that the Labware MiniHub has been loaded as shown in [Table 11](#).

**Table 11** Initial MiniHub configuration for LibraryPrep\_QXT\_ILM\_v1.0.rst

Vertical Shelf Position	Cassette 1	Cassette 2	Cassette 3	Cassette 4
Shelf 5 (Top)	Empty	Empty Nunc DeepWell plate	Empty	Empty
Shelf 4	Empty	Empty	Empty	Stop Solution source plate from <a href="#">step 13</a>
Shelf 3	Empty	Empty Eppendorf plate	Empty	Empty
Shelf 2	New tip box	Nuclease-free water reservoir from <a href="#">step 11</a>	AMPure XP beads in Nunc DeepWell plate from <a href="#">step 10</a>	Empty
Shelf 1 (Bottom)	Empty tip box	70% ethanol reservoir from <a href="#">step 12</a>	Empty	Empty tip box



## Step 2. Fragment and adaptor-tag the genomic DNA samples

20 Load the Bravo deck according to Table 12.

**Table 12** Initial Bravo deck configuration for LibraryPrep\_QXT\_ILM\_v1.0.rst


Location	Content
1	Empty waste reservoir (Axygen 96 Deep Well Plate, square wells)
4	Empty red insert
6	Library Prep Master Mix source plate (unsealed)
7	gDNA samples (5 $\mu$ L of 10 ng/ $\mu$ L DNA per well) in Eppendorf plate (unsealed)
9	Empty Eppendorf plate on red insert

### Run VWorks runset LibraryPrep\_QXT\_ILM\_v1.0.rst


21 On the SureSelect setup form, under **Select Protocol to Run**, select **LibraryPrep\_QXT\_ILM\_v1.0.rst**.

22 Select the number of columns of samples to be processed. Runs must include 1, 2, 3, 4, 6, or 12 columns.

23 Click **Display Initial Workstation Setup**.

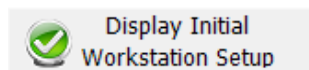

24 Verify that the NGS workstation has been set up as displayed in the Workstation Setup region of the form.

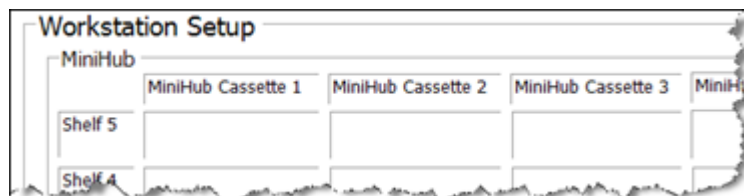


25 When verification is complete, click **Run Selected Protocol**.



### 3 Sample Preparation

#### Step 2. Fragment and adaptor-tag the genomic DNA samples

26 When ready to begin the run, click **OK** in the following window.

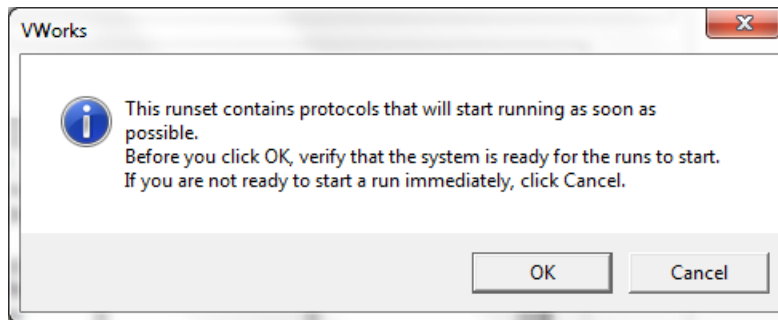

Running the LibraryPrep\_QXT\_ILM\_v1.0.rst runset takes approximately 1 hour. Once complete, the purified, adaptor-ligated DNA samples are located in the Eppendorf plate at position 7 of the Bravo deck.

## Step 3. Amplify adaptor-ligated libraries

In this step, the Agilent NGS Workstation completes the liquid handling steps for amplification and index tagging of the adaptor-ligated DNA samples using automation protocol WholeGenomePCR\_QXT\_ILM\_v1.0.pro. Afterward, you transfer the PCR plate to a thermal cycler for amplification.

This step uses the SureSelect<sup>QXT</sup> Reagent Kit components listed in [Table 13](#).

**Table 13** Reagents for whole genome library amplification

Kit Component	Storage Location	Where Used
Herculase II Fusion DNA Polymerase	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 45</a>
Herculase II 5× Reaction Buffer	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 45</a>
100 mM dNTP Mix (25 mM each dNTP)	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 45</a>
SureSelect QXT P7 and P5 dual indexing primers	SureSelect QXT Library Prep Kit Box 2, -20°C	<a href="#">page 46</a>
DMSO	Transferred to Room Temperature storage on <a href="#">page 34</a>	<a href="#">page 45</a>

### CAUTION

To avoid cross-contaminating libraries, set up PCR master mixes in a dedicated clean area or PCR hood with UV sterilization and positive air flow.

### Prepare the workstation

- 1 Clear the Labware MiniHub and BenchCel of all plates and tip boxes.
- 2 Turn on the ThermoCube, set to 0°C, at position 9 of the Bravo deck. Be sure that the chiller reservoir contains at least 300 mL of 25% ethanol.
- 3 Pre-set the temperature of Bravo deck positions 4 and 6 to 4°C using the Inheco Multi TEC control touchscreen, as described in [Setting the Temperature of Bravo Deck Heat Blocks](#). Bravo deck position 4 corresponds to CPAC 2, position 1 and deck position 6 corresponds to CPAC 2, position 2 on the Multi TEC control touchscreen.

### 3 Sample Preparation

#### Step 3. Amplify adaptor-ligated libraries

- 4 Load tip boxes for the run in the BenchCel Microplate Handling Workstation according to [Table 14](#).

**Table 14** Initial BenchCel configuration for WholeGenomePCR\_QXT\_ILM\_v1.0.pro

No. of Columns Processed	Rack 1	Rack 2	Rack 3	Rack 4
1	1 Tip box	Empty	Empty	Empty
2	1 Tip box	Empty	Empty	Empty
3	1 Tip box	Empty	Empty	Empty
4	1 Tip box	Empty	Empty	Empty
6	1 Tip box	Empty	Empty	Empty
12	1 Tip box	Empty	Empty	Empty

### Prepare the PCR master mix and master mix source plate

- 5 Prepare the appropriate volume of PCR Master Mix, according to [Table 15](#). Mix well using a vortex mixer and keep on ice.

**Table 15** Preparation of PCR Master Mix

SureSelect <sup>OXT</sup> Reagent	Volume for 1 Library	Volume for 1 Column	Volume for 2 Columns	Volume for 3 Columns	Volume for 4 Columns	Volume for 6 Columns	Volume for 12 Columns
Nuclease-free water	7.5 µL	95.6 µL	159.4 µL	223.1 µL	286.9 µL	414.4 µL	828.8 µL
Herculase II 5X Reaction Buffer	10.0 µL	127.5 µL	212.5 µL	297.5 µL	382.5 µL	552.5 µL	1105 µL
DMSO	2.5 µL	31.9 µL	53.1 µL	74.4 µL	95.6 µL	138.1 µL	276.3 µL
dNTP mix	0.5 µL	6.4 µL	10.6 µL	14.9 µL	19.1 µL	27.6 µL	55.3 µL
Herculase II Fusion DNA Polymerase	1.0 µL	12.8 µL	21.3 µL	29.8 µL	38.3 µL	55.3 µL	110.5 µL
<b>Total Volume</b>	<b>21.5 µL</b>	<b>274.1 µL</b>	<b>456.9 µL</b>	<b>639.6 µL</b>	<b>822.4 µL</b>	<b>1187.9 µL</b>	<b>2375.8 µL</b>

- 6 Using the same Nunc DeepWell master mix source plate that was used for the LibraryPrep\_QXT\_ILM\_v1.0.rst runset, add the volume of PCR master mix indicated in [Table 16](#) to all wells of column 2. Keep the source plate on ice until it is used on [page 47](#).

**Table 16** Preparation of the Master Mix Source Plate for WholeGenomePCR\_QXT\_ILM\_v1.0.pro

Master Mix Solution	Position on Source Plate	Volume of Master Mix added per Well of Nunc Deep Well Source Plate					
		1-Column Runs	2-Column Runs	3-Column Runs	4-Column Runs	6-Column Runs	12-Column Runs
PCR Master Mix	Column 2 (A2-H2)	31.6 µL	54.4 µL	77.3 µL	100.1 µL	145.8 µL	294.3 µL

### 3 Sample Preparation

#### Step 3. Amplify adaptor-ligated libraries

##### Assign and aliquot dual indexing primers

- 7 Determine the appropriate index assignments for each sample. See the [Reference](#) section for sequences of the index portion of the P7 and P5 indexing primers used to amplify the DNA libraries in this step. (See [Table 37](#) and [Table 38](#) for sequencing on HiSeq and MiSeq platforms or see [Table 37](#) and [Table 39](#) for sequencing on the NextSeq platform.)

Use the following guidelines for dual index assignments:

- Use a different indexing primer combination for each sample to be sequenced in the same lane.
  - All samples on the same row of the prepared DNA library plate must be assigned to the same P5 indexing primer (P5 i13 through P5 i20). This design results from the automation protocol configuration in which the P5 indexing primer is dispensed from a single source plate column to all columns of the indexing PCR plate. Each row of samples may be assigned to the same or different P5 primers, depending on run size and multiplexing requirements. (See [step 10](#), below, for P5 primer source plate setup details.)
  - The automation protocol configuration allows for any of the provided P7 indexing primers (P7 i1 through P7 i12) to be assigned to any sample position of the prepared DNA library plate. (See [step 8](#) and [step 9](#) below, for P7 primer source plate setup details.)
  - For sample multiplexing, Agilent recommends maximizing index diversity on both P7 and P5 primers as required for color balance. For example, when 8-plexing, use eight different P7 index primers with two P5 index primers. See [Table 40](#) on page 74 for additional details.
- 8 Dilute each P7 indexing primer (P7 i1 through P7 i12) to be used in the run according to [Table 17](#). The volumes below include the required excess.

**Table 17** Preparation of P7 indexing primer dilutions

Reagent	Volume to Index 1 Sample	Volume to Index 8 Samples
Nuclease-free water	3.0 µL	25.5 µL
SureSelect QXT P7 dual indexing primer (P7 i1 to P7 i12)	2.0 µL	17 µL
<b>Total Volume</b>	<b>5.0 µL</b>	<b>42.5 µL</b>

- 9 In a fresh PCR plate, aliquot 5  $\mu\text{L}$  of the appropriate P7 indexing primer dilution from [Table 17](#) to the intended sample indexing well position(s).

Keep the plate on ice.

- 10 Using the Nunc DeepWell master mix source plate containing PCR Master Mix from [step 6](#), add each P5 indexing primer (P5 i13 through P5 i20) to be used in the run to the appropriate well of column 2. Add the volume of P5 primer listed in [Table 18](#) to each well of column 2, according to the number of sample columns in the run. Each well of column 2 can contain the same or different P5 indexing primers.

The final configuration of the master mix source plate is shown in [Figure 5](#). Keep the source plate on ice.

**Table 18** Addition of P5 indexing primer to the DeepWell PCR source plate

Solution added to Source Plate	Position on Source Plate	Volume of Primer added per Well of Nunc Deep Well Source Plate					
		1-Column Runs	2-Column Runs	3-Column Runs	4-Column Runs	6-Column Runs	12-Column Runs
SureSelect QXT P5 dual indexing primer(s)*	Column 2 (A2-H2)	2.9 $\mu\text{L}$	5.1 $\mu\text{L}$	7.2 $\mu\text{L}$	9.3 $\mu\text{L}$	13.6 $\mu\text{L}$	27.4 $\mu\text{L}$


\* Each well of column 2 may contain the same or different P5 indexing primer. Typical 12-column runs include all eight of the provided SureSelect QXT P5 dual indexing primers (P5 i13 through P5 i20), resulting in a different P5 primer assignment to each row of the PCR indexing plate.

#### NOTE

If you are using a new DeepWell plate for the PCR source plate, leave column 1 empty and add the P5 Index(es) to column 2 of the new plate.

### 3 Sample Preparation

#### Step 3. Amplify adaptor-ligated libraries

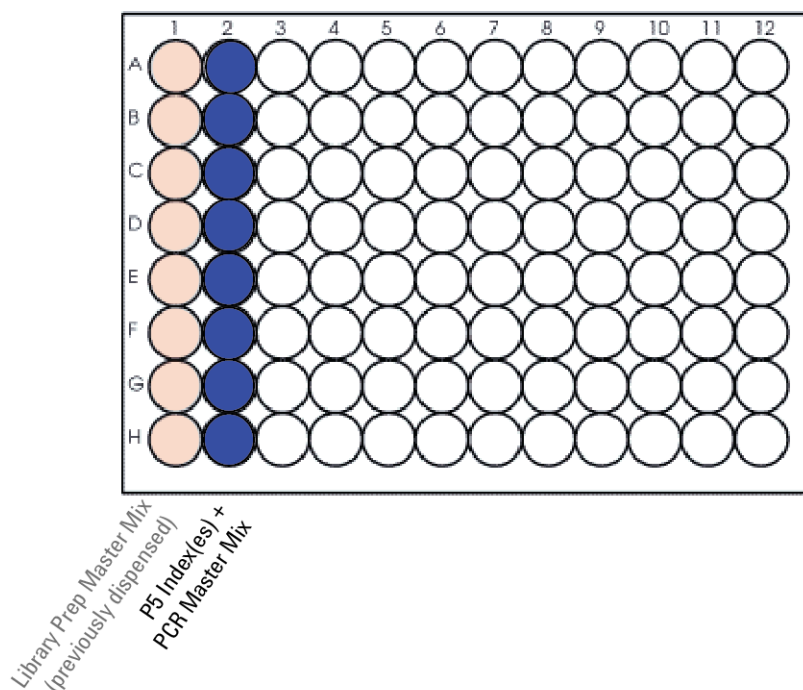

**Figure 5** Configuration of the master mix source plate for WholeGenomePCR\_QXT\_ILM\_v1.0.pro. Column 1 was used to dispense master mix during the previous protocol.

**11** Seal the master mix source plate using the PlateLoc Thermal Microplate Sealer, with sealing settings of 165°C and 1.0 sec.

**12** Centrifuge the plate for 30 seconds to drive the well contents off the walls and plate seal and to eliminate any bubbles.

#### NOTE

The presence of bubbles in source plate solutions may cause inaccurate volume transfer by the Bravo liquid handling platform. Ensure that the source plate is sealed and centrifuged prior to use in a run.



### Load the Agilent NGS Workstation

**13** Load the Labware MiniHub according to [Table 19](#).

**Table 19** Initial MiniHub configuration for WholeGenomePCR\_QXT\_ILM\_v1.0.pro

Vertical Shelf Position	Cassette 1	Cassette 2	Cassette 3	Cassette 4
Shelf 5 (Top)	Empty	Empty	Empty	Empty
Shelf 4	Empty	Empty	Empty	Empty
Shelf 3	Empty	Empty	Empty	Empty
Shelf 2	New tip box	Empty	Empty	Empty
Shelf 1 (Bottom)	Empty tip box	Empty	Empty	Empty tip box

**14** Load the Bravo deck according to [Table 20](#).

**Table 20** Initial Bravo deck configuration for WholeGenomePCR\_QXT\_ILM\_v1.0.pro

Location	Content
4	Adaptor-tagged DNA samples in Eppendorf plate
6	P7 indexing primers in PCR plate seated in red insert (PCR plate type must be specified on setup form under step 2)
9	Master mix plate containing P5 indexing primers and PCR Master Mix in Column 2 (unsealed)

### Run VWorks protocol WholeGenomePCR\_QXT\_ILM\_v1.0.pro

**15** On the SureSelect setup form, under **Select Protocol to Run**, select **WholeGenomePCR\_QXT\_ILM\_v1.0.pro**.


**16** Under **Select PCR plate labware for Thermal Cycling**, select the specific type of PCR plate used at position 6 of the Bravo deck.

**17** Select the number of columns of samples to be processed. Runs must include 1, 2, 3, 4, 6, or 12 columns.


### 3 Sample Preparation

#### Step 3. Amplify adaptor-ligated libraries


**18** Click **Display Initial Workstation Setup**.

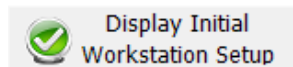

**19** Verify that the NGS workstation has been set up as displayed in the Workstation Setup region of the form.

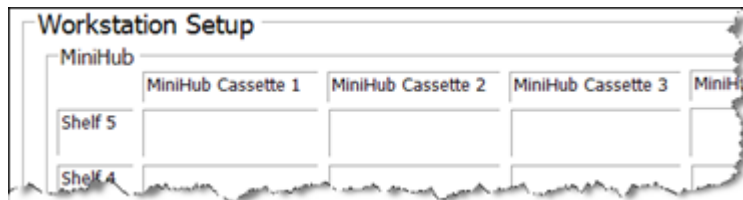


**20** When verification is complete, click **Run Selected Protocol**.

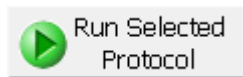

Running the WholeGenomePCR\_QXT\_ILM\_v1.0.pro protocol takes approximately 15 minutes. Once complete, the PCR-ready samples, containing prepped DNA, dual indexing primers, and PCR master mix, are located in the PCR plate at position 6 of the Bravo deck.

- 21** When you see the following prompt, remove the PCR plate from position 6 of the Bravo deck and seal the plate using the PlateLoc Thermal Microplate Sealer, with sealing settings of 165°C and 3.0 seconds.

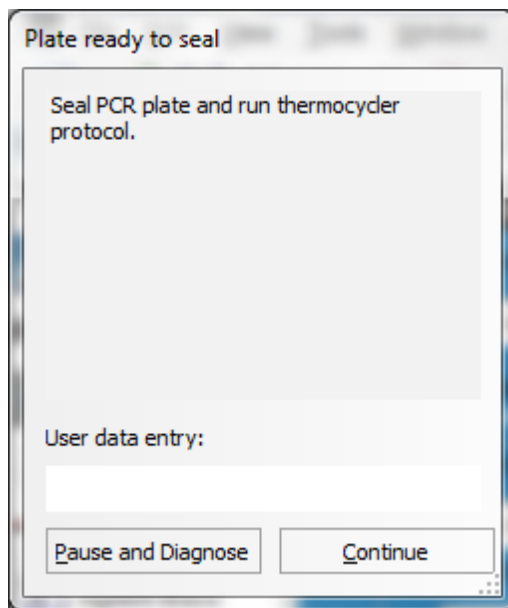

- 22** Centrifuge the plate for 30 seconds to drive the well contents off the walls and plate seal and to eliminate air bubbles.
- 23** Transfer the PCR plate to a thermal cycler and run the PCR amplification program shown in [Table 21](#).

### 3 Sample Preparation

#### Step 3. Amplify adaptor-ligated libraries

**Table 21** PCR cycling program

Segment	Number of Cycles	Temperature	Time
1	1	68°C	2 minutes
2	1	98°C	30 seconds
3	5	98°C	30 seconds
		56°C	30 seconds
		72°C	1 minute
4	1	4°C	Hold

## Step 4. Purify amplified DNA using AMPure XP beads

In this step, the Agilent NGS Workstation transfers AMPure XP beads to the indexed DNA sample plate and then collects and washes the bead-bound DNA.

### Prepare the workstation and reagents

- 1 Clear the Labware MiniHub and BenchCel of all plates and tip boxes.
- 2 Gently wipe down the Labware MiniHub, Bravo decks, and BenchCel with a DNA Away decontamination wipe.
- 3 Verify that the AMPure XP bead suspension is at room temperature. (If necessary, allow the bead solution to come to room temperature for at least 30 minutes.) *Do not freeze the beads at any time.*
- 4 Mix the bead suspension well so that the reagent appears homogeneous and consistent in color.
- 5 Turn on the ThermoCube, set to 4°C, at position 9 of the Bravo deck. Be sure that the chiller reservoir contains at least 300 mL of 25% ethanol.
- 6 Pre-set the temperature of Bravo deck position 4 to 45°C using the Inheco Multi TEC control touchscreen, as described in [Setting the Temperature of Bravo Deck Heat Blocks](#). Bravo deck position 4 corresponds to CPAC 2, position 1 on the Multi TEC control touchscreen.
- 7 Prepare a Nunc DeepWell source plate for the beads by adding 45 µL of homogeneous AMPure XP beads per well, for each well to be processed.
- 8 Prepare a Thermo Scientific reservoir containing 15 mL of nuclease-free water.
- 9 Prepare a separate Thermo Scientific reservoir containing 45 mL of freshly-prepared 70% ethanol.
- 10 Centrifuge the amplified DNA sample plate for 30 seconds to drive the well contents off the walls and plate seal and to remove air bubbles.

### 3 Sample Preparation

#### Step 4. Purify amplified DNA using AMPure XP beads

**11** Load the Labware MiniHub according to [Table 22](#), using the plate orientations shown in [Figure 2](#) on page 20.

**Table 22** Initial MiniHub configuration for AMPureXP\_QXT\_ILM\_v1.0.pro:Whole Genome PCR

Vertical Shelf Position	Cassette 1	Cassette 2	Cassette 3	Cassette 4
Shelf 5 (Top)	Empty Nunc DeepWell plate	Empty	Empty	Empty
Shelf 4	Empty	Empty	Empty	Empty
Shelf 3	Empty	Empty Eppendorf Plate	Empty	Empty
Shelf 2	Empty	Nuclease-free water reservoir from <a href="#">step 8</a>	AMPure XP beads in Nunc DeepWell plate from <a href="#">step 5</a>	Empty
Shelf 1 (Bottom)	Empty	70% ethanol reservoir from <a href="#">step 9</a>	Empty	Empty tip box

**12** Load the Bravo deck according to [Table 23](#).

**Table 23** Initial Bravo deck configuration for AMPureXP\_QXT\_ILM\_v1.0.pro:Whole Genome PCR

Location	Content
1	Empty waste reservoir (Axygen 96 Deep Well Plate, square wells)
9	Amplified DNA libraries in unsealed PCR plate seated in red insert (PCR plate type must be specified on setup form under step 2)

## Step 4. Purify amplified DNA using AMPure XP beads

**13** Load the BenchCel Microplate Handling Workstation according to Table 24.

**Table 24** Initial BenchCel configuration for AMPureXP\_QXT\_ILM\_v1.0.pro:Whole Genome PCR

No. of Columns Processed	Rack 1	Rack 2	Rack 3	Rack 4
1	1 Tip box	Empty	Empty	Empty
2	1 Tip box	Empty	Empty	Empty
3	2 Tip boxes	Empty	Empty	Empty
4	2 Tip boxes	Empty	Empty	Empty
6	3 Tip boxes	Empty	Empty	Empty
12	6 Tip boxes	Empty	Empty	Empty

**Run VWorks protocol *AMPureXP\_QXT\_ILM\_v1.0.pro:Whole Genome PCR***

**14** On the SureSelect setup form, under **Select Protocol to Run**, select **AMPureXP\_QXT\_ILM\_v1.0.pro:Whole Genome PCR**.


**NOTE**

AMPureXP purification protocols are used during multiple steps of the SureSelect automation workflow. Be sure to select the correct workflow step when initiating the automation protocol.

**15** Under **Select PCR plate labware for Thermal Cycling**, select the specific type of PCR plate containing the amplified libraries at position 9.

**16** Select the number of columns of samples to be processed. Runs must include 1, 2, 3, 4, 6, or 12 columns.


**17** Click **Display Initial Workstation Setup**.

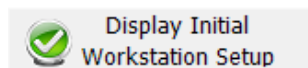

### 3 Sample Preparation

#### Step 4. Purify amplified DNA using AMPure XP beads

- 18 Verify that the NGS workstation has been set up as displayed in the Workstation Setup region of the form.

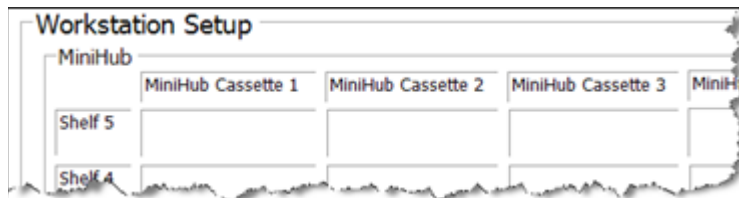

- 19 When verification is complete, click **Run Selected Protocol**.



The purification protocol takes approximately 45 minutes. When complete, the purified DNA samples are in the Eppendorf plate located on Bravo deck position 7.



## Step 5. Assess library DNA quantity and quality using the 2100 Bioanalyzer and High Sensitivity DNA Assay

Use the Bioanalyzer High Sensitivity DNA Assay to analyze the amplified indexed DNA. See the *High Sensitivity DNA Kit Guide* at [www.genomics.agilent.com](http://www.genomics.agilent.com) for more information on doing this step.

### NOTE

Do not use Agilent's DNA 1000 Assay to analyze the whole genome samples. The expected distribution of whole genome library fragment sizes is not compatible with the DNA 1000 Assay. See [Figure 6](#), below, for a sample electropherogram.

The presence of magnetic beads in the samples may adversely impact the Bioanalyzer results. If you suspect bead contamination in the samples, place the plate or strip tube on the magnetic rack before withdrawing samples for analysis.

- 1 Set up the 2100 Bioanalyzer as instructed in the reagent kit guide.
- 2 Prepare the chip, samples and ladder as instructed in the reagent kit guide, using 1  $\mu$ L of a 1:10 dilution of each sample for the analysis. Load the prepared chip into the 2100 Bioanalyzer and start the run within five minutes after preparation.
- 3 Analyze the results, using the guidelines below:
  - Typical whole genome library electropherograms show a broad distribution of DNA fragments. A sample electropherogram is shown in [Figure 6](#).
  - Check the Average Size [bp] of DNA fragments reported in the Bioanalyzer results. Sequencing data may be acquired from libraries with a broad range of average fragment sizes. The protocol has been optimized, however, to produce whole genome libraries with average DNA fragment sizes between approximately 600 and 1000 bp.

### NOTE


An average fragment size significantly less than 600 bp may indicate too little gDNA in the fragmentation reaction and may be associated with increased duplicates in the sequencing data. In contrast, libraries with an unusually large average fragment size may indicate too much gDNA in the fragmentation reaction and may require higher DNA concentrations for optimal cluster density in the sequencing reaction.

- 4 Measure the concentration of each library by integrating under the entire peak. For accurate quantification, make sure that the concentration falls within the linear range of the assay.

### 3 Sample Preparation

#### Step 5. Assess library DNA quantity and quality using the 2100 Bioanalyzer and High Sensitivity DNA Assay

**Stopping Point** If you do not continue to the next step, seal the plate and store at 4°C overnight or at -20°C for prolonged storage.

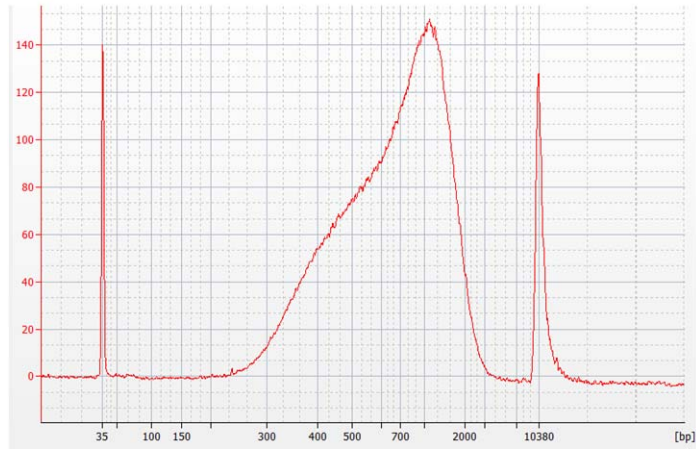

**Figure 6** Analysis of amplified library DNA using a High-Sensitivity DNA Assay.

## Step 6. Pool samples for multiplexed sequencing

The number of indexed libraries that may be multiplexed in a single sequencing lane is determined by the output specifications of the platform used, together with the amount of sequencing data required for your research design. Calculate the number of indexes that can be combined per lane, according to the capacity of your platform and the amount of sequencing data required per sample.

Guidelines for optimal low-level multiplexing of samples indexed using the SureSelect<sup>QXT</sup> dual indexes are provided on [page 74](#).

- 1 Combine the libraries such that each index-tagged sample is present in equimolar amounts in the pool. For each library, use the formula below to determine the amount of indexed sample to use.

$$\text{Volume of Index} = \frac{V(f) \times C(f)}{\# \times C(i)}$$

where  $V(f)$  is the final desired volume of the pool,

$C(f)$  is the desired final concentration of all the DNA in the pool

$\#$  is the number of indexes, and

$C(i)$  is the initial concentration of each indexed sample.

[Table 25](#) shows an example of the amount of 4 index-tagged samples (of different concentrations) and Low TE needed for a final volume of 20  $\mu\text{L}$  at 10 nM.

**Table 25** Example of indexed sample volume calculation for total volume of 20  $\mu\text{L}$

Component	V(f)	C(i)	C(f)	#	Volume to use ( $\mu\text{L}$ )
Sample 1	20 $\mu\text{L}$	20 nM	10 nM	4	2.5
Sample 2	20 $\mu\text{L}$	10 nM	10 nM	4	5
Sample 3	20 $\mu\text{L}$	17 nM	10 nM	4	2.9
Sample 4	20 $\mu\text{L}$	25 nM	10 nM	4	2
Low TE					7.6

- 2 Adjust the final volume of the pooled library to the desired final concentration.

### 3 Sample Preparation

#### Step 6. Pool samples for multiplexed sequencing

- If the final volume of the combined index-tagged samples is less than the desired final volume,  $V(f)$ , add Low TE to bring the volume to the desired level.
  - If the final volume of the combined index-tagged samples is greater than the final desired volume,  $V(f)$ , lyophilize and reconstitute to the desired volume.
- 3** If you store the library before sequencing, add Tween 20 to 0.1% v/v and store at  $-20^{\circ}\text{C}$  short term.

## Step 7. Prepare sequencing samples

The optimal seeding concentration for SureSelect<sup>QXT</sup> whole genome libraries is 14 to 20 pM on HiSeq or MiSeq instruments and 1.2 to 1.4 pM on the NextSeq platform. Seeding concentration and cluster density may also need to be optimized based on the DNA fragment size range for the library and on the desired output and data quality. Follow Illumina's recommendation for a PhiX control in a low-concentration spike-in for improved sequencing quality control.

### NOTE

The recommended seeding concentration of 14 to 20 pM is for samples quantified using the 2100 Bioanalyzer as described on [page 57](#). If you are using a different quantification method, such as QPCR, you may need to optimize the seeding concentration to achieve the optimal cluster density.

Proceed to cluster amplification using the appropriate Illumina Paired-End Cluster Generation Kit. See [Table 26](#) for kit configurations compatible with the recommended read length plus reads for the SureSelect<sup>QXT</sup> 8-bp dual indexes. To do this step, refer to the manufacturer's instructions, using the modifications described in "Using the SureSelect<sup>QXT</sup> Read Primers with Illumina's Paired-End Cluster Generation Kits" on page 62.

**Table 26** Illumina Kit Configuration Selection Guidelines

Platform	Run Type	Read Length*	SBS Kit Configuration	Chemistry
HiSeq 2500	Rapid Run	2 × 100 bp	200 Cycle Kit	v1 or v2
HiSeq 2500	High Output	2 × 100 bp	4 × 50 Cycle Kit <sup>†</sup>	v3
HiSeq 2500	High Output	2 × 125 bp	250 Cycle Kit	v4
HiSeq 2000	All Runs	2 × 100 bp	4 × 50 Cycle Kit <sup>†</sup>	v3
MiSeq	All Runs	2 × 150 bp	300 Cycle Kit	v2
MiSeq	All Runs	2 × 300 bp	600 Cycle Kit	v3
NextSeq 500	All Runs	2 × 150 bp	300 Cycle Kit	v2

\* If your application requires a different read length, verify that you have sufficient sequencing reagents to complete Reads 1 and 2 in addition to the dual 8-bp index reads.

† A single 200-cycle kit does not include enough reagents to complete Reads 1 and 2 in addition to the dual 8-bp index reads in this format. If preferred, the additional reads may be supported by using one 200-cycle kit plus one 50-cycle kit.

### 3 Sample Preparation

#### Step 7. Prepare sequencing samples

#### Using the SureSelect<sup>QXT</sup> Read Primers with Illumina's Paired-End Cluster Generation Kits

To sequence the SureSelect<sup>QXT</sup> libraries on Illumina sequencing platforms, you need to use the following custom sequencing primers provided in SureSelect QXT Library Prep Kit Box 2:

- SureSelect QXT Read Primer 1
- SureSelect QXT Read Primer 2
- SureSelect QXT Index Read Primer
- SureSelect QXT Index 2 Read Primer NSQ (NextSeq platform only)

These SureSelect<sup>QXT</sup> custom sequencing primers are provided at 100  $\mu\text{M}$  and must be diluted 1:200 in the corresponding Illumina primer solution, using the platform-specific instructions below:

**For the HiSeq platform**, combine the primers as shown in [Table 27](#) or [Table 28](#) on [page 63](#).

**For the MiSeq platform**, combine the primers as shown in [Table 29](#) on [page 63](#).

**For the NextSeq platform**, combine the primers as shown in [Table 30](#) or [Table 31](#) on [page 64](#).

#### NOTE

It is important to combine the primers precisely in the indicated ratios. Be sure to use measured volumes of each solution; do not use volumes reported on vial labels when preparing the mixtures. Vortex each mixture vigorously to ensure homogeneity for proper detection of the indexes using the custom read primers.

**Table 27 HiSeq2000 and HiSeq 2500 High Output custom sequencing primer preparation**

Sequencing Read	Volume of SureSelect QXT Primer	Volume of Illumina TruSeq Primer	Total Volume
Read 1	5 µL SureSelect QXT Read Primer 1 (brown cap)	995 µL HP6 or HP10	1 mL*
Index	15 µL SureSelect QXT Index Read Primer (clear cap)	2985 µL HP8 or HP12	3 mL
Read 2	15 µL SureSelect QXT Read Primer 2 (black cap)	2985 µL HP7 or HP11	3 mL

\* Aliquot the mixture as directed for HP6 or HP10 in Illumina's cluster generation protocol.

**Table 28 HiSeq 2500 Rapid Mode custom sequencing primer preparation**

Sequencing Read	Volume of SureSelect QXT Primer	Volume of Illumina TruSeq Primer	Total Volume
Read 1	8.8 µL SureSelect QXT Read Primer 1 (brown cap)	1741.2 µL HP10	1.75 mL*
Index	8.8 µL SureSelect QXT Index Read Primer (clear cap)	1741.2 µL HP12	1.75 mL
Read 2	8.8 µL SureSelect QXT Read Primer 2 (black cap)	1741.2 µL HP11	1.75 mL

\* Aliquot the mixture as directed for HP10 in Illumina's cluster generation protocol.

**Table 29 MiSeq platform custom sequencing primer preparation**

Sequencing Read	Volume of SureSelect <sup>QXT</sup> Primer	Volume of Illumina TruSeq Primer	Total Volume	Final Cartridge Position
Read 1	3 µl SureSelect QXT Read Primer 1 (brown cap)	597 µl HP10 (well 12)	0.6 ml	well 18
Index	3 µl SureSelect QXT Index Read Primer (clear cap)	597 µl HP12 (well 13)	0.6 ml	well 19
Read 2	3 µl SureSelect QXT Read Primer 2 (black cap)	597 µl HP11 (well 14)	0.6 ml	well 20

### 3 Sample Preparation

#### Step 7. Prepare sequencing samples

**Table 30** NextSeq 500/550 High-Output v2 Kit custom sequencing primer preparation

Sequencing Read	Volume of SureSelect <sup>QXT</sup> Primer	Volume of Illumina Primer	Total Volume	Final Cartridge Position
Read 1	3.9 µl SureSelect QXT Read Primer 1 (brown cap)	1296.1 µl BP10 (from well 20)	1.3 ml	well 7
Read 2	4.2 µl SureSelect QXT Read Primer 2 (black cap)	1395.8 µl BP11 (from well 21)	1.4 ml	well 8
Index + Index 2	6 µl SureSelect QXT Index Read Primer (clear cap) + 6 µl SureSelect QXT Index 2 Read Primer NSQ (purple cap)	1988 µl BP14 (from well 22)	2 ml	well 9

**Table 31** NextSeq 500/550 Mid-Output v2 Kit custom sequencing primer preparation

Sequencing Read	Volume of SureSelect <sup>QXT</sup> Primer	Volume of Illumina Primer	Total Volume	Final Cartridge Position
Read 1	2.7 µl SureSelect QXT Read Primer 1 (brown cap)	897.3 µl BP10 (from well 20)	0.9 ml	well 7
Read 2	3.3 µl SureSelect QXT Read Primer 2 (black cap)	1096.7 µl BP11 (from well 21)	1.1 ml	well 8
Index + Index 2	4.8 µl SureSelect QXT Index Read Primer (clear cap) + 4.8 µl SureSelect QXT Index 2 Read Primer NSQ (purple cap)	1590.4 µl BP14 (from well 22)	1.6 ml	well 9



## Step 8. Set up the sequencing run and trim adaptors from the reads

Refer to Illumina protocols to set up custom sequencing primer runs, using the additional guidelines outlined below.

For SureSelect<sup>QXT</sup> dual index sequence information, see tables on [page 72](#) through [page 73](#).

Before aligning reads to the reference genome, SureSelect<sup>QXT</sup> adaptor sequences must be trimmed from the reads. You can use SureCall, Agilent's NGS data analysis software, to perform adaptor trimming, alignment of reads and variant calling of sequencing data generated from either the HiSeq or the MiSeq platform. To download SureCall free-of-charge and for additional information, including tutorials on this software, visit the [SureCall page at www.genomics.agilent.com](http://www.genomics.agilent.com).

SureCall is compatible with FASTQ files generated by both the HiSeq and MiSeq platforms. To use SureCall to analyze SureSelect<sup>QXT</sup>-generated data, you first need to define an analysis workflow. This analysis workflow identifies the libraries as SureSelect<sup>QXT</sup> libraries and enables automated adaptor trimming. The trimmed FASTQ files can then be used for alignment to generate BAMs for downstream analysis.

To create the analysis workflow, refer to [Figure 7 on page 66](#). Upon starting SureCall, click the **Analysis Workflow** tab. Choose the appropriate analysis type (single sample, paired, or trio analysis), and then click the **Import Unaligned Files** button. Within the *Select Unaligned Sample Files* window, specify your read 1 and read 2 files using the **Add** buttons. Using the menus near the bottom of the screen, select **Default SureSelect QXT Method** from the *Analysis Method* menu, choose the appropriate design description from the *Design* menu, and select **Illumina** from the *Platform* menu. Once done, refer to the SureCall guide for next steps on alignment and variant calling.

If using another pipeline for alignment and downstream analysis, refer to the platform-specific guidelines starting on [page 66](#).

### 3 Sample Preparation

#### Step 8. Set up the sequencing run and trim adaptors from the reads

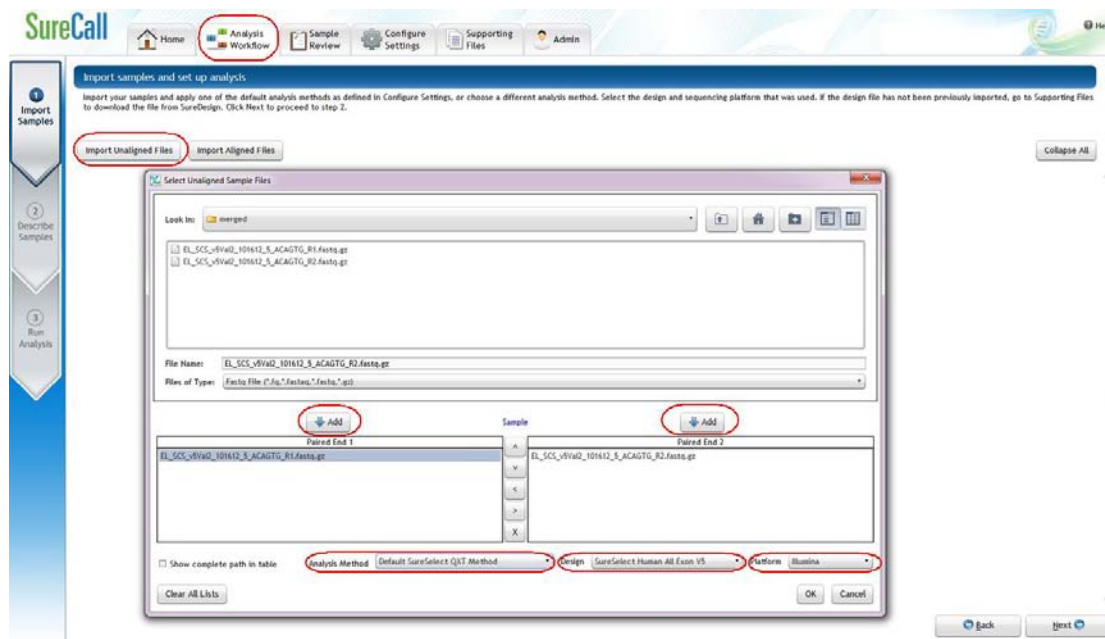


Figure 7 Analysis workflow setup in SureCall.

#### MiSeq platform sequencing run setup and adaptor trimming guidelines

Use the Illumina Experiment Manager (IEM) software to generate a custom primer Sample Sheet.

Set up the run to include adaptor trimming using the IEM Sample Sheet Wizard. When prompted by the wizard, select the *Use Adaptor Trimming* option, and specify **CTGTCCTTGATCACA** as the adaptor sequence. This enables the MiSeq Reporter software to identify the adaptor sequence and trim the adaptor from reads.

#### HiSeq or NextSeq 500 platform sequencing run setup and adaptor trimming guidelines

Set up sequencing runs using the *Custom* setting. Since custom primers are spiked into the standard sequencing primer tubes, no additional specialized settings are required to accommodate the use of custom primers in the run.

## Step 8. Set up the sequencing run and trim adaptors from the reads

Use the *Cycles* settings shown in Table 32. Cycle number settings can be specified on the *Run Configuration* screen of the instrument control software interface after choosing *Custom* from the index type selection buttons.

**Table 32** Run Configuration screen Cycle Number settings

Run Segment	Cycle Number
Read 1	100
Index 1 (i7)	8
Index 2 (i5)	8
Read 2	100

After the sequencing run is complete, generate demultiplexed FASTQ data following Illumina's instructions and then trim adaptor sequences from the reads using Agilent's Read Trimmer tool. This tool takes in data in FASTQ format and removes the adaptor sequence from the ends of the sequencing reads, generating trimmed FASTQ data as output. To download the Read Trimmer tool free-of-charge and for additional information on this resource, visit [www.agilent.com/genomics](http://www.agilent.com/genomics).

### **3 Sample Preparation**

**Step 8. Set up the sequencing run and trim adaptors from the reads**



## 4 Reference

Kit Contents	70
Nucleotide Sequences of SureSelect <sup>OXT</sup> Dual Indexes	72
Guidelines for Multiplexing with Dual-Indexed Samples	74

This chapter contains reference information, including component kit contents and reference information for use during the downstream sample sequencing steps.



## Kit Contents

SureSelect<sup>QXT</sup> Library Prep Kits contain the following component kits:

**Table 33** SureSelect<sup>QXT</sup> Library Prep Kit Contents

Component Kits	Storage Condition	HiSeq or MiSeq (ILM) 96 Samples	NextSeq (NSQ) 96 Samples
SureSelect QXT Library Prep Kit, ILM, Box 1	Room Temperature	5500-0119	5500-0119
SureSelect QXT Library Prep Kit Box 2	-20°C	5500-0121	5500-0127

The contents of each of the component kits listed in [Table 33](#) are described in [Table 34](#) through [Table 36](#) below.

**Table 34** SureSelect QXT Library Prep, ILM, Box 1 Content

Kit Component	96 Reactions
SureSelect QXT Stop Solution	bottle

**Table 35** SureSelect QXT Library Prep, ILM, Box 2 Content

Kit Component	HiSeq or MiSeq (ILM) 96 Reactions
SureSelect QXT Buffer	bottle
SureSelect QXT Enzyme Mix ILM	tube with orange cap
Herculase II Fusion DNA Polymerase	tube with red cap
Herculase II 5× Reaction Buffer	tube with clear cap
100 mM dNTP Mix (25 mM each dNTP)	tube with green cap
DMSO	tube with green cap
SureSelect QXT Read Primer 1	tube with amber cap
SureSelect QXT Read Primer 2	tube with black cap
SureSelect QXT Index Read Primer	tube with clear cap
SureSelect QXT P7 dual indexing primers	P7 i1 through P7 i12 provided in 12 tubes with yellow caps (one tube per primer)
SureSelect QXT P5 dual indexing primers	P5 i13 through P5 i20 provided in 8 tubes with blue caps (one tube per primer)

**Table 36** SureSelect QXT Library Prep, NSQ, Box 2 Content

Kit Component	NextSeq (NSQ) 96 Reactions
SureSelect QXT Buffer	bottle
SureSelect QXT Enzyme Mix ILM	tube with orange cap
Herculase II Fusion DNA Polymerase	tube with red cap
Herculase II 5× Reaction Buffer	tube with clear cap
100 mM dNTP Mix (25 mM each dNTP)	tube with green cap
DMSO	tube with green cap
SureSelect QXT Read Primer 1	tube with amber cap
SureSelect QXT Read Primer 2	tube with black cap
SureSelect QXT Index Read Primer	tube with clear cap
SureSelect QXT Index 2 Read Primer NSQ	tube with purple cap
SureSelect QXT P7 dual indexing primers	P7 i1 through P7 i12 provided in 12 tubes with yellow caps (one tube per primer)
SureSelect QXT P5 dual indexing primers	P5 i13 through P5 i20 provided in 8 tubes with blue caps (one tube per primer)

## Nucleotide Sequences of SureSelect<sup>QXT</sup> Dual Indexes

The nucleotide sequence of each SureSelect<sup>QXT</sup> index is provided in the tables below.

Note that some index number assignments of the SureSelect<sup>QXT</sup> P5 and P7 indexes differ from the index numbers assignments used by Illumina for indexes of similar or identical sequence.

Each index is 8 bases in length. Refer to Illumina's sequencing run setup instructions for sequencing libraries using 8-base indexes.

**Table 37** SureSelect<sup>QXT</sup> P7 Indexes 1 to 12

Index Number	Sequence
P7 Index 1 (P7 i1)	TAAGGCGA
P7 Index 2 (P7 i2)	CGTACTAG
P7 Index 3 (P7 i3)	AGGCAGAA
P7 Index 4 (P7 i4)	TCCTGAGC
P7 Index 5 (P7 i5)	GTAGAGGA
P7 Index 6 (P7 i6)	TAGGCATG
P7 Index 7 (P7 i7)	CTCTCTAC
P7 Index 8 (P7 i8)	CAGAGAGG
P7 Index 9 (P7 i9)	GCTACGCT
P7 Index 10 (P7 i10)	CGAGGCTG
P7 Index 11 (P7 i11)	AAGAGGCA
P7 Index 12 (P7 i12)	GGACTCCT



**Table 38** SureSelect<sup>OXT</sup> P5 Indexes 13 to 20 for HiSeq platform, MiSeq platform, or NextSeq platform runs through BaseSpace

Index Number	Sequence
P5 Index 13 (P5 i13)	TAGATCGC
P5 Index 14 (P5 i14)	CTCTCTAT
P5 Index 15 (P5 i15)	TATCCTCT
P5 Index 16 (P5 i16)	AGAGTAGA
P5 Index 17 (P5 i17)	GTAAGGAG
P5 Index 18 (P5 i18)	ACTGCATA
P5 Index 19 (P5 i19)	AAGGAGTA
P5 Index 20 (P5 i20)	CTAAGCCT

**Table 39** SureSelect<sup>OXT</sup> P5 Indexes 13 to 20 for NextSeq platform\*

Index Number	Sequence
P5 Index 13 (P5 i13)	GCGATCTA
P5 Index 14 (P5 i14)	ATAGAGAG
P5 Index 15 (P5 i15)	AGAGGATA
P5 Index 16 (P5 i16)	TCTACTCT
P5 Index 17 (P5 i17)	CTCCTTAC
P5 Index 18 (P5 i18)	TATGCAGT
P5 Index 19 (P5 i19)	TACTCCTT
P5 Index 20 (P5 i20)	AGGCTTAG

\* When doing NextSeq runs through BaseSpace, use the reverse complement sequences provided in [Table 38](#).

## 4 Reference

### Guidelines for Multiplexing with Dual-Indexed Samples

## Guidelines for Multiplexing with Dual-Indexed Samples

Agilent recommends following the dual index sample pooling guidelines shown in Table 40. These are designed to maintain color balance at each cycle of the index reads on both ends. They also provide flexibility of demultiplexing as single or dual indexed samples in low-plexity experiments. One-base mismatches should also be allowed during demultiplexing in order to maximize sequencing output per sample.

**Table 40** Dual index sample pooling guidelines for 96 Reaction Kits

Plexity of Sample Pool	Recommended SureSelect <sup>QXT</sup> P7 Indexes	Recommended SureSelect <sup>QXT</sup> P5 Indexes
1-plex	Any P7 index i1 to i11	Any P5 index (i13 to i20)
2-plex	P7 i1 and P7 i2 OR P7 i2 and P7 i4	P5 i13 and P5 i14 OR P5 i15 and P5 i16 OR P5 i17 and P5 i18
3-plex	P7 i1, P7 i2 and P7 i4 OR P7 i3, P7 i4 and P7 i6 OR P7 i5, P7 i7 and P7 i8	P5 i13 and P5 i14 OR P5 i15 and P5 i16 OR P5 i17 and P5 i18 (as needed)
4-plex	P7 i1, P7 i2, P7 i3* and P7 i4 OR P7 i3, P7 i4, P7 i5* and P7 i6 OR P7 i5, P7 i6*, P7 i7 and P7 i8	P5 i13 and P5 i14 OR P5 i15 and P5 i16 OR P5 i17 and P5 i18 (as needed)
5-plex	P7 i1, P7 i2, P7 i3*, P7 i4 and P7 i5* OR P7 i3, P7 i4, P7 i5*, P7 i6 and p7 i7* OR P7 i5, P7 i6*, P7 i7, P7 i8 and p7 i9*	P5 i13 and P5 i14 OR P5 i15 and P5 i16 OR P5 i17 and P5 i18 (as needed)
6- to 11-plex	Any combination of P7 indexes i1 to i11 using each index only once	P5 i13 and P5 i14 OR P5 i15 and P5 i16 OR P5 i17 and P5 i18 (as needed)
12- to 88-plex	Any combination of P7 indexes i1 to i11 (as needed)	P5 i13 and P5 i14 and any third P5 index OR P5 i15 and P5 i16 and any third P5 index OR P5 i17 and P5 i18 and any third P5 index (as needed)
89- to 96-plex	All twelve P7 indexes (i1 to i12)	Any P5 indexes (i13 to i20, as needed)

\* The indicated indexes may be substituted with another index within range of P7 i1 to P7 i11, as long as the substitute index differs from all others used in the sample pool.

[www.agilent.com](http://www.agilent.com)

## In This Book

This guide contains information to run the SureSelect<sup>QXT</sup> Automated Library Prep protocol.

© Agilent Technologies, Inc. 2014, 2015

Version C0, November 2015



p/n G9682-90010



**Agilent Technologies**