



***Taq* Extender PCR Additive**

Instruction Manual

Catalog #600148 (1000 U)

Revision C.0

For Research Use Only. Not for use in diagnostic procedures.

600148-12



Agilent Technologies

LIMITED PRODUCT WARRANTY

This warranty limits our liability to replacement of this product. No other warranties of any kind, express or implied, including without limitation, implied warranties of merchantability or fitness for a particular purpose, are provided by Agilent. Agilent shall have no liability for any direct, indirect, consequential, or incidental damages arising out of the use, the results of use, or the inability to use this product.

ORDERING INFORMATION AND TECHNICAL SERVICES

Email

techservices@agilent.com

World Wide Web

www.genomics.agilent.com

Telephone

Location	Telephone
United States and Canada	800 227 9770
Austria	01 25125 6800
Benelux	02 404 92 22
Denmark	45 70 13 00 30
Finland	010 802 220
France	0810 446 446
Germany	0800 603 1000
Italy	800 012575
Netherlands	020 547 2600
Spain	901 11 68 90
Sweden	08 506 4 8960
Switzerland	0848 8035 60
UK/Ireland	0845 712 5292
All Other Countries	Please visit www.agilent.com/genomics/contactus

Taq Extender PCR Additive

CONTENTS

Materials Provided	1
Storage Conditions	1
Additional Materials Required	1
Notices To Purchaser	1
Introduction	2
Protocol	2
Troubleshooting	3
Preparation of Reagents	3
References	3
Endnotes	3

Taq Extender PCR Additive

MATERIALS PROVIDED

Materials provided	Quantity
Taq Extender PCR additive (5 U/ μ l)	200 μ l
Taq Extender 10 \times reaction buffer ^a	1 ml

^a See Preparation of Reagents.

STORAGE CONDITIONS

All components: -20°C

ADDITIONAL MATERIALS REQUIRED

Taq DNA polymerase

NOTICES TO PURCHASER

Limited Label License: The purchase of this product conveys to the purchaser only the limited, non-transferable right to use the purchased quantity of the product for the purchaser's own research by the purchaser only under patent claims outside the U.S. corresponding to the expired U.S. Patent No. 5,436,149 (claims 6-16). No rights are granted to the purchaser to sell, modify for resale or otherwise transfer this product, either alone or as a component of another product, to any third party. Takara Bio reserves all other rights, and this product may not be used in any manner other than as provided herein. For information on obtaining a license to use this product for purposes other than research, please contact Takara Bio Inc., Seta 3-4-1, Otsu, Shiga 520-2193, Japan (Fax +81-77-543-9254).

NOTICE TO PURCHASER: LIMITED LICENSE

Use of this product is covered by one or more of the following US patents and corresponding patent claims outside the US: 6,258,569, 6,171,785, 6,127,155, 6,030,787, 5,994,056, 5,876,930, 5,804,375, 5,789,224, 5,773,258 (claims 1 and 6 only), 5,723,591, 5,677,152 (claims 1 to 23 only), 5,618,711, 5,538,848, and claims outside the US corresponding to expired US Patent No. 5,079,352. The purchase of this product includes a limited, non-transferable immunity from suit under the foregoing patent claims for using only this amount of product for the purchaser's own internal research. No right under any other patent claim and no right to perform commercial services of any kind, including without limitation reporting the results of purchaser's activities for a fee or other commercial consideration, is conveyed expressly, by implication, or by estoppel. This product is for research use only. Diagnostic uses under Roche patents require a separate license from Roche. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

INTRODUCTION

Taq Extender PCR additive is a polymerase enhancer that improves the reliability and yield of conventional polymerase chain reaction (PCR) amplifications. This additive increases the efficiency at which *Taq* DNA polymerase performs extension reactions on specific DNA segments in each cycle of PCR, thus resulting in a greater percentage of the extension reactions reaching completion. In addition, *Taq* Extender PCR additive improves the PCR amplification of difficult templates and increases the reliability and yield of many PCR targets up to 10 kb in length.¹ The easy-to-use *Taq* Extender PCR additive is simply added to amplification reactions in an amount equal to that of *Taq* DNA polymerase, the standard *Taq* 10× reaction buffer is replaced with an optimized *Taq* Extender 10× reaction buffer and cycling is performed using standard PCR conditions.

PROTOCOL

1. Prepare the amplification reaction using a standard PCR protocol and implement the modifications to the standard protocol as described in step 2. Determine the standard reaction parameters, such as the deoxynucleotide (dNTP) concentration, the amount of *Taq* DNA polymerase required and the cycling conditions (see step 3).

2. Modify the standard PCR protocol as outlined below.

Notes *Modifications of primer length or cycling temperatures are not necessary.*

*Use of thin-wall microcentrifuge tubes is highly recommended, but not critical, to improve the heat transfer, which will further enhance the efficacy of the *Taq* Extender PCR additive.*

- a. Substitute the standard *Taq* 10× reaction buffer with the *Taq* Extender 10× reaction buffer.
 - b. Add an equal number of units of *Taq* Extender PCR additive and *Taq* DNA polymerase to the amplification reaction [i.e., add 1 µl of *Taq* Extender PCR additive (5 U/µl) to an amplification reaction requiring 1 µl of *Taq* DNA polymerase (5 U/µl)]. Use a minimum of 1 U each of *Taq* DNA polymerase and *Taq* Extender PCR additive per kilobase to be amplified.
3. Perform the amplification reaction using standard cycling conditions, allowing at least 30–60 seconds of extension time for each kilobase to be amplified.²

TROUBLESHOOTING

Observation	Solution(s)
Low yield	Minor lot-to-lot variations in the concentration of <i>Taq</i> DNA polymerases from various manufacturers may affect PCR product yields and may necessitate the use of greater or lesser amounts of <i>Taq</i> Extender PCR additive per amplification reaction in order to achieve optimal PCR results
	Choice of the <i>Taq</i> Extender 10× reaction buffer may also affect PCR product yields. If the <i>Taq</i> Extender 10× reaction buffer provided with this kit does not appear to be optimized for your particular amplification system, further buffer optimization may be required
	Insufficient amounts of <i>Taq</i> DNA polymerase or inadequate extension times may also contribute to poor PCR product yields. Optimal PCR product yields are obtained by using a minimum of 1 U of <i>Taq</i> DNA polymerase per kilobase to be amplified and by allowing at least 30–60 seconds of extension time for each kilobase to be amplified ²
Artificial smears	PCR smearing may be due primarily to minor contaminants of the <i>Taq</i> DNA polymerase. Two possible solutions are to use a different lot number of the enzyme or to reduce the extension time

PREPARATION OF REAGENTS

Taq Extender 10× Reaction Buffer

- 200 mM Tris-HCl (pH 8.8)
- 100 mM KCl
- 100 mM (NH₄)₂SO₄
- 20 mM MgSO₄
- 1% Triton® X-100
- 1 mg/ml nuclease-free bovine serum albumin (BSA)

REFERENCES

1. Nielson, K., Schoettlin, W., Bauer, J. C. and Mathur, E. (1994) *Strategies* 7(2):27.
2. Innis, M. A., Myambo, K. B., Gelfand, D. H. and Brow, M. A. (1988) *Proc Natl Acad Sci U S A* 85(24):9436-40.

ENDNOTES

Triton® is a registered trademark of Rohm and Haas Co.

MSDS INFORMATION

Material Safety Data Sheets (MSDSs) are provided online at <http://www.genomics.agilent.com>. MSDS documents are not included with product shipments.